

User-Developed Protocol:

Isolation of genomic and viral DNA from lymphocytes using the QIAGEN® Genomic-tip

This procedure has been adapted by customers from the Genomic-tip Protocol, and is for use with the QIAGEN® Genomic-tip. **It has not been thoroughly tested or optimized by QIAGEN.**

Please be sure to read the *QIAGEN Genomic DNA Handbook* and the detailed Cell Cultures Protocol carefully before beginning this procedure.

Procedure

1. Centrifuge lymphocytes to form a pellet. Discard the supernatant.
2. Resuspend pellet in Buffer C1 as described in the Sample Preparation and Lysis Protocol for Cell Cultures in the *QIAGEN Genomic DNA Handbook*.
3. Centrifuge the lysed cells at 4°C for 15 min at 1300 x g. Save the nuclear pellet and the supernatant (S1).

Viral DNA in cytosolic fraction

- 3a. To supernatant S1, add RNase A to a final concentration of 10–20 µg/ml. Add 25 µl QIAGEN Protease or Proteinase K per ml of supernatant.
- 4a. Incubate for 1 h at 50°C.
- 5a. Add NaCl to a final concentration of 750 mM and adjust the pH to 7.0.
- 6a. Load onto an equilibrated QIAGEN Genomic-tip and follow the wash and elution steps in the Genomic-tip Protocol for Isolation of Genomic DNA from Blood, Cultured Cells, Tissue, Yeast, or Bacteria in the *QIAGEN Genomic DNA Handbook*.

Note: QIAGEN Genomic-tip 20/G should be used for the viral DNA purification steps. If the expected yield of viral DNA is unknown or less than 1 µg, a carrier, for example glycogen, should be added to the eluate before the isopropanol precipitation.

Viral DNA in nuclear fraction — integrated

- 3b. Treat nuclear pellet from Step 2 with Buffers P1, P2, and P3 and centrifuge as described in the QIAGEN Midi and Maxi Kits Protocols in the *QIAGEN Plasmid Purification Handbook*. Save the pellet (which contains the genomic DNA) and the supernatant (S2).
- 4b. Resuspend the pellet in Buffer G2, add Proteinase K or QIAGEN Protease, and incubate according to the instructions in Step 7 of the Sample Preparation and Lysis Protocol for Cell Cultures of the *QIAGEN Genomic DNA Handbook*.

**User-developed
protocol**

- 5b. Precipitate the DNA with 0.7 volumes of isopropanol and resuspend DNA in TE. Add Buffer QBT to a final volume of 1 ml for a Genomic-tip 20/G, 10 ml for a Genomic-tip 100/G, or 25 ml for a Genomic-tip 500/G. Load the sample onto the corresponding equilibrated Genomic-tip and follow the wash and elution steps in the *QIAGEN Genomic DNA Handbook* (Step 3 and Step 4 of the Protocol for Isolation of Genomic DNA from Blood, Cultured Cells, Tissue, Yeast, or Bacteria),

Viral DNA in nuclear fraction — not integrated

- 3c. Load S2 from Step 3b onto the equilibrated Genomic-tip and follow the wash and elution steps in the *QIAGEN Genomic DNA Handbook*.

Note: QIAGEN Genomic-tip 20/G should be used for the viral DNA purification steps. If the expected yield of viral DNA is unknown or less than 1 µg, a carrier, for example glycogen, should be added to the eluate before the isopropanol precipitation.

QIAGEN handbooks can be requested from QIAGEN Technical Service or your local QIAGEN distributor. Selected handbooks can be downloaded from www.qiagen.com/literature/handbooks/default.asp. Material safety data sheets (MSDS) for any QIAGEN product can be downloaded from www.qiagen.com/ts/msds.asp.

Trademark: QIAGEN® (QIAGEN).

© 2001 QIAGEN, all rights reserved.