

**User-developed
protocol**

User-Developed Protocol:

Isolation of bacterial DNA from soil using the QIAamp[®] DNA Stool Mini Kit and QIAamp DNA Blood Midi Kit

This procedure has been adapted by customers from the QIAamp[®] DNA Stool Mini Kit Protocols and is for use with the QIAamp DNA Stool Mini Kit and QIAamp DNA Blood Midi Kit. **It has not been thoroughly tested and optimized by QIAGEN.**

As starting material, 5 g soil was mixed with different amounts of *Bacillus subtilis* cells. Sensitivity was 5×10^3 cells/5g soil.

Please be sure to read the *QIAamp DNA Stool Mini Kit Handbook* carefully before beginning this procedure.

Procedure

DNA isolation

1. Weigh up to 5 g soil in a 50 ml BD Falcon™ tube.
2. Add 2–5 ml distilled water to the tube, and mix for 5 min on a shaker.
3. Incubate for 10 min at 95°C.
4. Centrifuge at 3000 rpm for 5 min. Transfer the supernatant to a new tube.
5. Add 7 volumes of Buffer ASL to the supernatant, and mix well
6. Add 1 InhibitEX™ tablet to the tube and incubate for 1 min at room temperature (15–25°C) on a shaker.
7. Centrifuge sample at 5000 x g for 5 min. Transfer the supernatant into a new tube.
8. Add 1 volume of Buffer AL to the supernatant, and mix well.
9. Add 1 volume of ethanol (96–100%).
10. Place a QIAamp Midi Spin Column on the QIAvac 24 vacuum manifold.
11. Apply the sample lysate onto the QIAamp Midi Spin Column. Apply maximum vacuum.
12. Wash the column once with 1 ml Buffer AW1.
13. Wash the column once with 1 ml Buffer AW2.
14. Place the QIAamp Midi Spin Column in a 15 ml tube (provided), and centrifuge at 5000 rpm for 15 min to dry the membrane.

**User-developed
protocol**

15. Place the QIAamp Midi Spin Column in a clean 15 ml tube. To elute the DNA, add 300 µl Buffer AE, and centrifuge at 5000 rpm for 5 min.
16. Reload the eluate onto the membrane of the QIAamp Midi Spin Column, and centrifuge at 5000 rpm for 5 min.

Amplification of a 528 bp *atpase* gene fragment from *B. subtilis*

Primer	Sequence 5'-3'
UEB1	GTGTGATTGTTTTATTGATTGC
UEB2	GTACCGACAAGACCGAGAGC

PCR mix

25 µl 10x HotStarTaq™ Master Mix
1 µl primer UEB1, 10 µM
1 µl primer UEB2, 10 µM
22 µl water
1 µl DNA (eluate)

Amplification conditions

95°C for 15 min	1x
94°C for 1 min; 52°C for 1 min 45 s; 72°C for 3 min	50x
72°C for 10 min	1x

For up-to-date licensing information and product-specific disclaimers, see the respective QIAGEN kit handbook or user manual. QIAGEN kit handbooks and user manuals are available at www.qiagen.com or can be requested from QIAGEN Technical Services or your local distributor.

QIAGEN handbooks can be requested from QIAGEN Technical Service or your local QIAGEN distributor. Selected handbooks can be downloaded from www.qiagen.com/literature/handbooks/default.aspx. Material safety data sheets (MSDS) for any QIAGEN product can be downloaded from www.qiagen.com/ts/msds.asp.

Trademarks: QIAGEN®, QIAamp®, QIAvac, HotStarTaq™; InhibitEX™ (QIAGEN); BD Falcon™ (Becton, Dickinson and Company).

QA28 © 2002–2010 QIAGEN, all rights reserved.