

QIASymphony[®] — Pure Performance




Sample & Assay Technologies



Relax and let QIAGEN® automate your complete workflow

QIAGEN now provides everything that you need to standardize and fully automate your routine sample preparation and analysis workflow. With our dedicated range of QIAAsymphony Kits, the [QIAAsymphony SP](#) enables sample preparation of DNA, RNA, bacterial and viral nucleic acids, and 6xHis-tagged protein from a wide range of starting materials. The [QIAAsymphony AS](#) extends the capabilities of the [QIAAsymphony SP](#) by integrating automated PCR assay setup which, in combination with the [Rotor-Gene® Q](#) and QIAGEN [real-time and end-point PCR kits](#), enables you to complete your automated PCR workflow and maximize your efficiency — from sample to result.

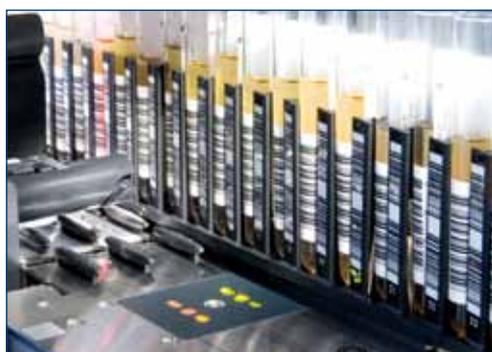

► The QIAAsymphony provides:

- Standardized sample preparation and PCR assay setup
- In-built touchscreen and prefilled reagent cartridges for ease of use
- Liquid-level detection, active cooling, and UV lamp for decontamination
- Applications for certified QIAGEN and in-house assays
- Service for laboratory-developed tests*

* For customization service for in-house assays, please inquire.




Cutting-edge system for standardized sample preparation and PCR setup

The [QIASymphony](#) system is ideal for laboratories that perform sample processing and PCR analysis on a day-to-day basis. With predefined protocols and automatable QIAGEN kits for sample preparation and PCR setup, the [QIASymphony](#) system will optimize your workflow, increasing the productivity in your laboratory. The [QIASymphony SP](#) is already the first choice for many laboratories, enabling success in virus and bacterial analysis, gene expression analysis, and in forensics and genomics applications. The high level of standardization that the [QIASymphony](#) system achieves allows you to directly compare your results with laboratories around the world. QIAGEN offers tailored IQ/OQ products to support laboratory certification for full instrument qualification.*

Modular system for maximum convenience and process safety

Laboratories already operating a [QIASymphony SP](#) can easily upgrade their system onsite with a [QIASymphony AS](#) module. The [QIASymphony AS](#) directly interfaces with the [QIASymphony SP](#), delivering fast, highly reproducible PCR setup. To reduce manual handling steps and minimize the risk of sample contamination, samples processed on the [QIASymphony SP](#) can be transferred automatically to the [QIASymphony AS](#) (integrated operation). For extra flexibility, the [QIASymphony SP and AS](#) can also be operated independently.† Samples that have been processed earlier in the day, or that have been processed on a different [QIASymphony SP](#), can be manually transferred to the [QIASymphony AS](#).

QIASymphony workflows

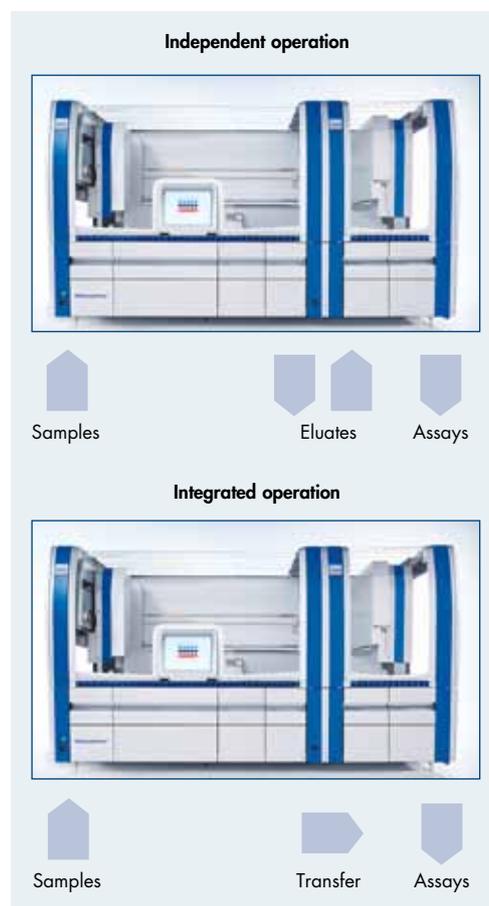

* For a full list of QIASymphony service agreements, please inquire. † The QIASymphony SP and AS remain physically connected during independent operation.

Ethernet adapter

For connecting the QIASymphony system to local networks

Touchscreen

For intuitive operation of the complete system

USB port

For bi-directional data exchange

Sample drawer

Accommodates 1–96 sample tubes plus internal controls in a tube carrier, or up to 4 microtiter plates in a plate carrier

Reagent and consumables drawer

Supports consumables and 2 reagent cartridges

Waste drawer

Stores solid and liquid waste separately

Eluate drawer


Provides a choice of elution formats, including cooling and transfer of eluate racks to the QIASymphony AS from the rear position


UV lamp

Decontaminates the worktable and minimizes sample cross-contamination

Robotic arm

4-channel with tip guards

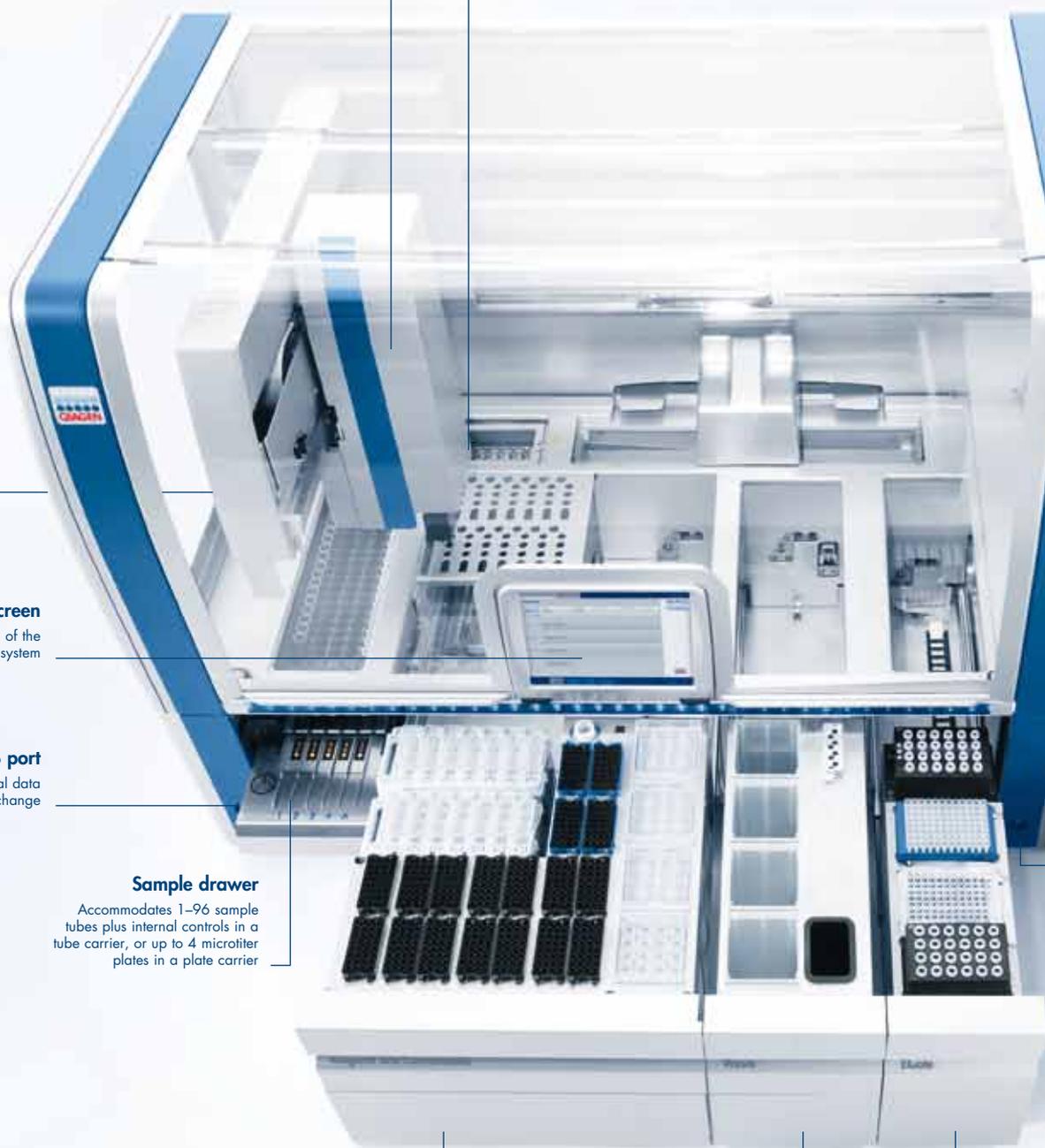

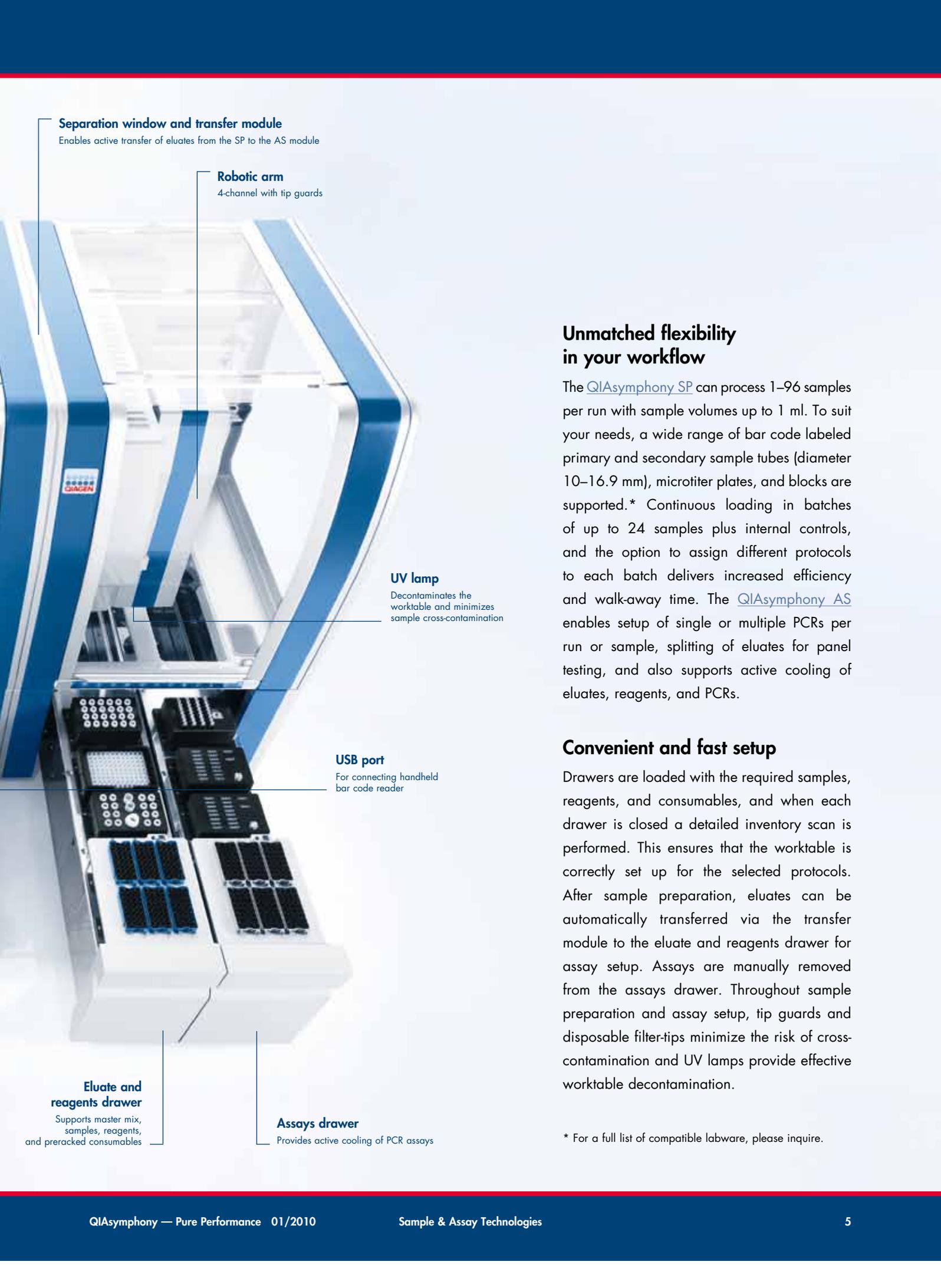

Separation window and transfer module

Enables active transfer of eluates from the SP to the AS module

Robotic arm

4-channel with tip guards

UV lamp

Decontaminates the worktable and minimizes sample cross-contamination

USB port

For connecting handheld bar code reader

Eluate and reagents drawer

Supports master mix, samples, reagents, and preracked consumables

Assays drawer

Provides active cooling of PCR assays

Unmatched flexibility in your workflow

The [QIASymphony SP](#) can process 1–96 samples per run with sample volumes up to 1 ml. To suit your needs, a wide range of bar code labeled primary and secondary sample tubes (diameter 10–16.9 mm), microtiter plates, and blocks are supported.* Continuous loading in batches of up to 24 samples plus internal controls, and the option to assign different protocols to each batch delivers increased efficiency and walk-away time. The [QIASymphony AS](#) enables setup of single or multiple PCRs per run or sample, splitting of eluates for panel testing, and also supports active cooling of eluates, reagents, and PCRs.

Convenient and fast setup

Drawers are loaded with the required samples, reagents, and consumables, and when each drawer is closed a detailed inventory scan is performed. This ensures that the worktable is correctly set up for the selected protocols. After sample preparation, eluates can be automatically transferred via the transfer module to the eluate and reagents drawer for assay setup. Assays are manually removed from the assays drawer. Throughout sample preparation and assay setup, tip guards and disposable filter-tips minimize the risk of cross-contamination and UV lamps provide effective worktable decontamination.

* For a full list of compatible labware, please inquire.



Sample information after sample loading

Easy-to-use software

[QIASymphony](#) operating software guides the user through setting up a sample preparation or assay setup run in a step-by-step fashion, with options to select and adapt standardized protocols. Required volumes of buffers, reagents, master mix, and controls are automatically calculated and the user is informed of the number and type of consumables to be loaded. Transfer of sample loading information from the [QIASymphony SP](#) to the [QIASymphony AS](#) is convenient and reduces the number of steps required for assay definition.



Assay definition of 2 PCR assays

Standardized data exchange permits LIMS compatibility

The [QIASymphony](#) system provides standardized communication interfaces. Work lists and rack files can be imported to the [QIASymphony](#) system, enabling automatic run definition before samples are received in the laboratory. After assay setup, sample lists can be exported to selected real-time PCR cyclers, including the [Rotor-Gene Q](#). This bi-directional data exchange allows integration of [QIASymphony](#) systems into laboratory information management systems (LIMS).



Reagent loading information for the [QIASymphony AS](#)

Seamless sample traceability

Bar code reading of samples and reagents enables full sample tracking throughout the entire purification and assay setup procedure. The [QIASymphony](#) system provides chronological electronic documentation of results, including information about sample IDs, their positions on the worktable, and the lot number and expiry date of sample preparation and PCR reagents. Multi-level user management provides a record of which user was responsible for sample processing. Electronic result files can be automatically printed via a networked printer, or can be downloaded using a USB stick for archiving purposes.

► Unrivaled application range on the QIAasymphony SP

QIAasymphony Kits — backed by QIAGEN Quality® — increase laboratory efficiency since just a few kits in combination with optimized protocols cover an extensive range of starting materials and applications (Table 1). Sample volumes of up to 1 ml, and up to 50 mg of tissue or 10^7 cells can be processed, providing increased nucleic acid yields for gene expression and genomics applications and higher sensitivity in virology and microbiology applications.

The QIAasymphony SP delivers pure performance in virus and bacteria applications and makes no compromise when purifying DNA from blood, RNA from tissues and cells, or His-tagged proteins from *E. coli* or eukaryotic cells. For up-to-date information about QIAasymphony applications, visit www.qiagen.com/goto/QIAasymphony.



Innovative reagent cartridges for ultra-efficient purification

QIAasymphony Kits are safe and exceptionally easy to use. Ready-to-run, bar code labeled reagent cartridges are prefilled with all reagents required for the purification procedure, including accessory enzymes. Worktable setup is rapid and saves you valuable time. Simply place up to 2 reagent cartridges in the QIAasymphony drawer. The reagent cartridges can be either from the same kit or different kits, enabling different purification procedures to be performed within the same run of 96 samples. Reagent cartridges are automatically opened by the instrument and reagent volumes only for the selected number of samples are used, giving you complete cost control.



Table 1. Dedicated QIAasymphony Kits for a wide range of starting materials

| QIAasymphony Kit | Starting material and amount |
|------------------------------------|--|
| ► QIAasymphony Virus/Bacteria | Purification of viral nucleic acids from up to 1000 μ l plasma, serum, or CSF. Purification of viral nucleic acids and bacterial DNA from up to 800 μ l respiratory samples, urine, stool, transport media, or air-dried swabs |
| ► QIAasymphony DNA | Purification of DNA from up to 1000 μ l human whole blood, 400 μ l buffy coat, 50 mg tissue, or 1×10^7 cells |
| ► QIAasymphony DNA Investigator | Purification of reference and casework samples for human identity testing and forensics |
| ► QIAasymphony AXpH DNA | Purification of DNA from liquid-based cytology samples using AXpH technology |
| ► QIAasymphony RNA | Purification of total RNA from up to 50 mg tissue or 1×10^7 cells |
| ► QIAasymphony PAXgene® Blood RNA* | Purification of cellular RNA from 2.5 ml human whole blood collected in PAXgene Blood RNA Tubes |
| ► QIAasymphony Ni-NTA Native | Purification of 6xHis-tagged proteins from up to 5 ml <i>E. coli</i> or up to 2.5×10^7 eukaryotic cells |
| ► QIAasymphony Ni-NTA Denaturing | Purification of 6xHis-tagged proteins from up to 5 ml <i>E. coli</i> or up to 2.5×10^7 eukaryotic cells |

* Up to 48 samples can be processed per run.

Table 3. Real-time PCR detection of HIV RNA from plasma

| Input titer (IU/ml) | Number of replicates | Number of positives | Percentage of positives |
|---------------------|----------------------|---------------------|-------------------------|
| 600 | 12 | 12 | 100 |
| 200 | 14 | 14 | 100 |
| 100 | 14 | 14 | 100 |
| 75 | 14 | 14 | 100 |
| 50 | 14 | 13 | 92.9 |
| 25 | 14 | 10 | 71.4 |
| 12.5 | 14 | 5 | 35.7 |
| 6 | 14 | 6 | 42.9 |
| 3 | 14 | 1 | 7.1 |
| 0 | 8 | 0 | 0 |

Serial dilutions from a negative plasma pool spiked with WHO HIV-1 RNA 2nd International Standard NIBSC Code: 97/650 were processed in replicates of 8–14 per titer using the [QIASymphony SP](#) with the Virus Cellfree 1000 protocol (60 µl elution volume) and the [QIASymphony Virus/Bacteria Midi Kit](#). HIV RNA was detected using an HI Virus-1 RT-PCR assay.

Table 4. Real-time PCR detection of CMV DNA from plasma

| Input titer (Copies/ml) | Number of replicates | Number of positives | Percentage of positives |
|-------------------------|----------------------|---------------------|-------------------------|
| 1000 | 24 | 24 | 100 |
| 316 | 24 | 24 | 100 |
| 100 | 24 | 24 | 100 |
| 32 | 24 | 23 | 95.8 |
| 10 | 24 | 15 | 62.5 |
| 3.2 | 24 | 4 | 16.7 |
| 1 | 24 | 2 | 8.3 |
| 0.32 | 24 | 1 | 4.2 |
| 0 | 9 | 0 | 0 |

Serial dilutions from a negative plasma pool spiked with CMV material with known titer were processed in replicates of 24 or 9 per titer using the [QIASymphony SP](#) with the Virus Cellfree 1000 protocol (60 µl elution volume) and the [QIASymphony Virus/Bacteria Midi Kit](#). CMV DNA was detected using a CMV PCR assay.

Purify viral nucleic acids or bacterial DNA

High-quality viral DNA and RNA, and bacterial nucleic acids can be purified from a wide range of sample types ([Tables 2–4](#)) using [QIASymphony Virus/Bacteria Kits](#) on the [QIASymphony SP](#). Optimized protocols deliver high yields of pure nucleic acids, and flexible elution volumes support your downstream assays. The [QIASymphony SP](#) is provided with standardized sample preparation protocols that define sample transfer, purification, and then elution. For purification of difficult samples or inactivation of samples, additional lysis procedures can be combined with the purification protocols.

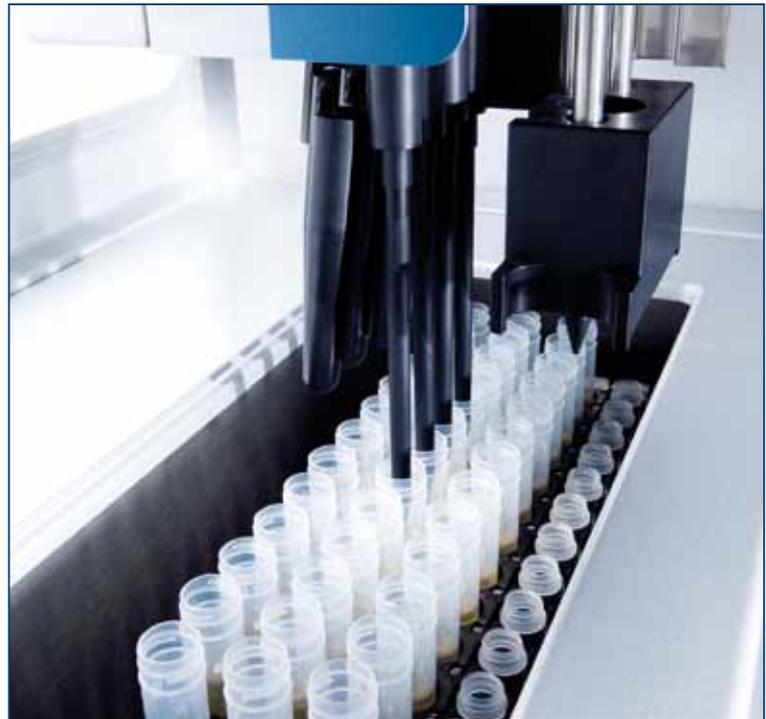


Table 2. Successful purification of a wide range of respiratory viruses from different sample types using the [QIASymphony SP](#)

| Respiratory Virus | Sample type |
|------------------------|---|
| Adenovirus | aspirates, swab |
| HSV1 | swab |
| Influenza A/B | aspirates, BAL, sputum, swab, pharyngeal lavage |
| hMPV | nasal secretion |
| Parainfluenzavirus 1–4 | aspirates, BAL, nasal secretion, swab |
| Coronavirus | aspirates, nasal secretion, sputum, swab |
| Rhinovirus | nasal secretion, tracheal and pharyngeal lavage |
| Enterovirus | nasal secretion, pharyngeal lavage |
| RSV | aspirates, sputum, swab, nasal secretion, pharyngeal lavage |
| Bocavirus | nasal secretion, tracheal and pharyngeal lavage |

Unmatched yield and purity

[QIAasymphony DNA Kits](#) enable automated purification of DNA from a broad range of sample types with sample volumes up to 1 ml. Genomic DNA can be purified from human and animal whole blood ([Figure 1](#)), buffy coat, human and animal tissues (up to 50 mg), cultured cells (up to 1×10^7), and bacterial cultures (up to 4×10^9 cells) ([Table 5](#)). The [QIAasymphony DNA Mini Kit](#) is also suited for purification of viral DNA from human whole blood ([Table 6](#)). There is a dedicated [QIAasymphony DNA Investigator Kit](#) for processing forensic reference and casework samples.

Optimized protocols provide high yields of pure, inhibitor-free DNA. Flexible elution volumes allow DNA concentration to be easily adapted to your needs. Purified DNA is ready for use in even the most demanding downstream applications, such as multiplex ligation-dependent probe amplification (MLPA), array comparative genomic hybridization (aCGH), SNP analyses, Southern hybridizations, and STR profiling ([Figure 2](#)).

Table 5. Different DNA yields obtained from a range of sample types

| Sample type | Sample amount or volume | Elution volume (µl) | Typical DNA yield (µg) |
|--------------|-------------------------|---------------------|------------------------|
| Whole blood | 200 µl | 200 | 4–12 |
| | 400 µl | 400 | 8–24 |
| | 1000 µl | 500 | 15–45 |
| Buffy coat | 200 µl | 200 | 12–40 |
| | 400 µl | 400 | 24–72 |
| Spleen | 25 mg | 200 | 40–80 |
| Liver | 25 mg | 200 | 25–50 |
| Muscle | 50 mg | 200 | 5–15 |
| Lung | 25 mg | 200 | 10–25 |
| Kidney | 25 mg | 200 | 15–30 |
| Rat tail | 50 mg | 200 | 20–40 |
| Jurkat cells | 1×10^7 cells | 200 | 60–80 |

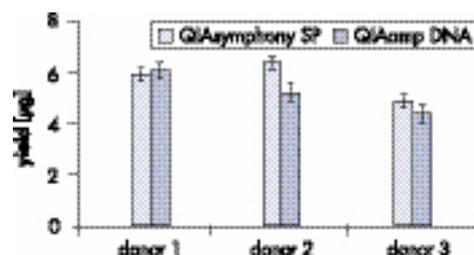
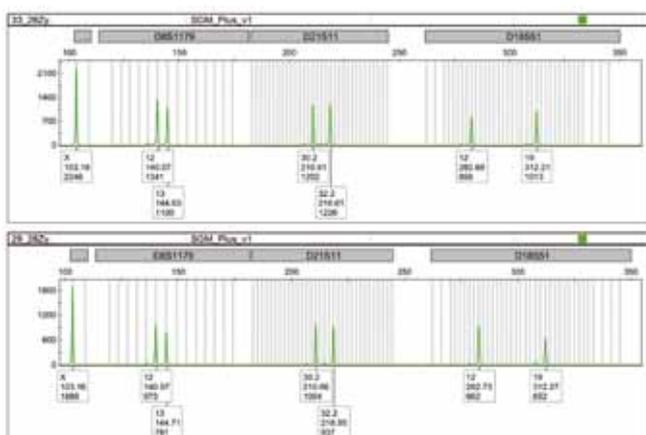


Figure 1. Equivalent performance to QIAamp. DNA was purified from 200 µl human whole blood. Replicates of 6, from 3 blood donors, were processed on the [QIAasymphony SP](#) using the [QIAasymphony DNA Mini Kit](#) and the [QIAamp DNA Blood Mini Kit](#).

Table 6. Real-time PCR detection of CMV DNA in human whole blood

| Input titer (Copies/ml) | Number of replicates | Number of positives | Percentage of positives |
|-------------------------|----------------------|---------------------|-------------------------|
| 250 | 33 | 33 | 100 |
| 200 | 33 | 33 | 100 |
| 150 | 39 | 38 | 97 |
| 100 | 45 | 40 | 89 |
| 75 | 12 | 8 | 67 |
| 50 | 45 | 27 | 60 |
| 25 | 26 | 6 | 23 |
| 10 | 24 | 5 | 21 |
| 0 | 15 | 0 | 0 |

Dilutions of CMV stock (Acrometrix) were processed in replicates of 12–45 using the [QIAasymphony SP](#) with the Virus Blood 200 protocol (60 µl elution volume) and the [QIAasymphony DNA Mini Kit](#). CMV DNA was detected using a CMV PCR assay.

Figure 2. STR profiling using the QIAasymphony SP and the EZ1 Advanced. DNA from saliva (10 µl) was purified on the [QIAasymphony SP](#) using the Casework 200 µl protocol and the [QIAasymphony DNA Investigator Kit](#), or on the [EZ1 Advanced](#) using the Trace protocol and the [EZ1 DNA Investigator Kit](#). DNA was eluted in 50 µl of Buffer ATE. DNA (1 ng) was analyzed by short tandem repeat (STR) profiling in 12.5 µl reactions. Data are shown for ameloginin, D8S1179, D21S11, and D18S51 loci.

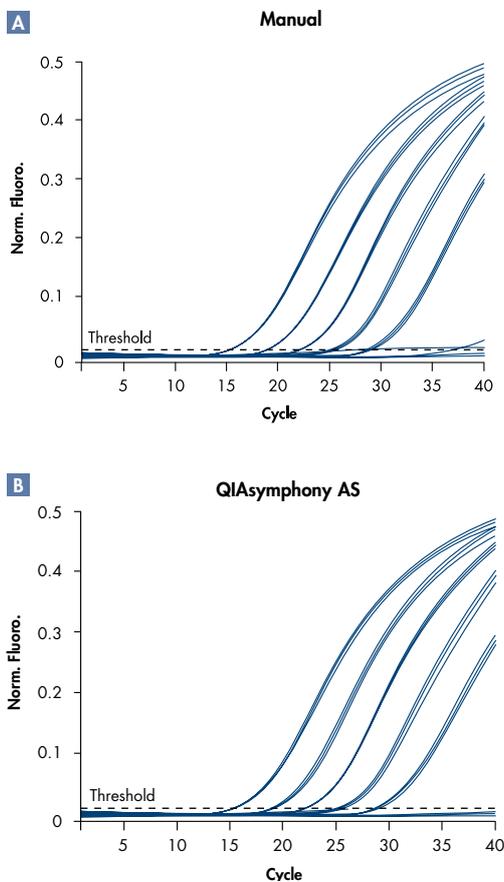


Figure 3. Highly reproducible assay setup. A series dilution of leukocyte cDNA (100 ng to 0.01 ng) was performed **A** manually and **B** using the [QIASymphony AS](#). GAPDH was amplified from these dilutions, using the [Rotor-Gene Q](#) and the [QuantiFast® Probe PCR Kit](#). Reactions were set up in replicates of 4 in Strip Tubes.

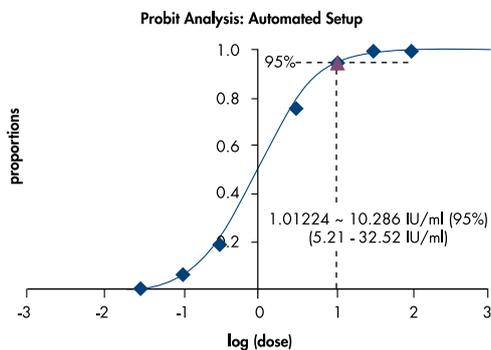


Figure 4. Automated assay setup is as sensitive as manual assay setup. Half-log dilutions of WHO HBV standard (n=8) were performed in replicates of 8. Assay setup was performed manually and using the [QIASymphony AS](#). Amplification was performed using the [Rotor-Gene Q](#) and a validated, in-house real-time PCR assay. Automated and manual assay setup were equally sensitive. Manual setup: 1,04698 ~ 11,142 IU/ml, (6,039–33,22 IU/ml)

Transfer your purified samples directly to the [QIASymphony AS](#)

The [QIASymphony AS](#) can simplify your routine work processes — even setting up complex PCR assays is fast and easy. Manual pipetting steps that are prone to human error are eliminated, so the [QIASymphony AS](#) delivers more accurate pipetting on a day-to-day basis than even highly experienced laboratory staff ([Figure 3 and 4](#)). Pipetting of master mix and sample transfer for 96 samples takes less than 25 minutes, leaving you and your staff with more valuable time for other tasks.

The [QIASymphony AS](#) is provided with predefined protocols, specifically designed for use with QIAGEN's expanding portfolio of [real-time, end-point, and *artus*® PCR kits](#) ([Table 7](#)). Standardized pipetting schemes can be adapted as required by QIAGEN customization services. A choice of output adapters enables use of different real-time PCR cyclers (e.g., [Rotor-Gene Q](#), 96-well cyclers, 32-capillary cyclers) for detection. You can now relax and let QIAGEN's automated solutions do the hard work for you — prepare your samples on the [QIASymphony SP](#) and then directly transfer your eluates to the [QIASymphony AS](#) for assay setup. To use the [Rotor-Gene Q](#) for detection, simply set up your PCRs in Strip Tubes or Rotor-Discs. Your workflow has never been easier!

Table 7. Automate proven QIAGEN PCR kits with the [QIASymphony AS](#)*

| Application | Method | Kit | Description |
|----------------------------|--------------------------|-----------------------------------|---|
| ▶ Gene expression analysis | Real-time RT-PCR | ▶ Rotor-Gene Kits | Rapid and reliable RT-PCR on the Rotor-Gene Q |
| | Real-time RT-PCR | ▶ QuantiFast Kits | Fast, sensitive RT-PCR on any real-time cycler |
| | Real-time RT-PCR | ▶ QuantiTect® Kits | Sensitive RT-PCR on any real-time cycler |
| | End-point RT-PCR | ▶ QIAGEN OneStep RT-PCR Kit | Sensitive RT-PCR, with simple setup, using any RNA template |
| ▶ Virus detection | Real-time PCR and RT-PCR | ▶ QuantiTect Kits | Highly sensitive detection of multiple viral DNA and RNA targets in the same tube |
| | End-point RT-PCR | ▶ QIAGEN OneStep RT-PCR Kit | Highly sensitive detection of viral RNA |
| ▶ Pathogen detection | Real-time PCR and RT-PCR | ▶ <i>cador</i> ™ veterinary | Ready-to-use, validated real-time PCR assays to detect veterinary pathogens |
| | Real-time PCR and RT-PCR | <i>artus</i> PCR and RT-PCR Kits† | Ready-to-use real-time PCR assays for pathogen detection |

* For a full list of QIAGEN PCR kits, including kits for genotyping and other specialized applications, visit www.qiagen.com/goto/PCR. For availability of [QIASymphony AS](#) protocols, please inquire. † *cador* veterinary PCR and RT-PCR Kits are not available in the US or Canada. ‡ *artus* PCR and RT-PCR Kits are not available in the US.

Ordering Information

| Product | Contents | Cat. no. |
|--|---|----------|
| ▶ QIAAsymphony SP | QIAAsymphony sample prep module, 1-year warranty on parts and labor | 9001297 |
| ▶ QIAAsymphony AS | QIAAsymphony assay setup module, 1-year warranty on parts and labor | 9001301 |
| ▶ Rotor-Gene Q 5plex HRM | Real-time PCR cyclers and High Resolution Melt analyzer with 5 channels | 9001580 |
| ▶ QIAAsymphony Cabinet SP/AS | For correct positioning of the QIAAsymphony SP/AS instruments | 9020246 |
| ▶ Starter Pack, QIAAsymphony AS | Pack includes consumables required for operating the QIAAsymphony AS | 997199 |
| ▶ Filter-Tips, 50 µl, Qsym AS (1024) | Disposable Filter-Tips; racked; (8 x 128) | 997120 |
| ▶ QIAAsymphony Virus/Bacteria Mini Kit (192) | For up to 192 preps of 200 µl each: Includes 2 reagent cartridges and enzyme racks and accessories | 931036 |
| ▶ QIAAsymphony Virus/Bacteria Midi Kit (96) | For up to 96 preps of 1000 µl each: Includes 2 reagent cartridges and enzyme racks and accessories | 931055 |
| ▶ QIAAsymphony DNA Mini Kit (192) | For 192 preps of 200 µl each: Includes 2 reagent cartridges and enzyme racks and accessories | 931236 |
| ▶ QIAAsymphony DNA Midi Kit (96) | For 96 preps of 1000 µl each: Includes 2 reagent cartridges and enzyme racks and accessories | 931255 |
| ▶ QIAAsymphony AXpH DNA Kit (192) | For 192 preps: Includes 2 reagent cartridges and enzyme racks and accessories | 937156 |
| ▶ QIAAsymphony DNA Investigator Kit | For 192 preps from casework and reference samples: Includes 2 reagent cartridges and enzyme racks and accessories | 931436 |
| ▶ QIAAsymphony RNA Kit (192) | For 192 preps: Includes 2 reagent cartridges and enzyme racks and accessories | 931636 |
| ▶ QIAAsymphony PAXgene Blood RNA Kit (96) | For 962 preps: Includes 2 reagent cartridges and enzyme racks and accessories | 762535 |

For up-to-date licensing information and product-specific disclaimers, see the respective QIAGEN kit handbook or user manual. QIAGEN kit handbooks and user manuals are available at www.qiagen.com or can be requested from QIAGEN Technical Services or your local distributor.

Visit www.qiagen.com/goto/QIAAsymphony to find out more!

Trademarks: QIAGEN®, QIAGEN Quality®, QIAamp®, QIAAsymphony®, artus®, EZ1®, QuantiFast®, QuantiTect® (QIAGEN Group); Rotor-Disc™, Rotor-Gene® (Corbett Research); PAXgene® (PreAnalytiX GmbH).

The QIAGEN silver logo is exclusively licensed to Corbett Research.

NOTICE TO PURCHASER:

The purchase of this product (Rotor-Gene Q, Rotor-Disc) includes a limited, non-transferable license to certain patents (see details below) surrounding rapid polymerase chain reaction (PCR) methods and instrumentation, the use of SYBR® Green I in PCR reactions, melting curve analysis, analysis methods of DNA melting data, specifically high resolution melting (HRM) and others.

The purchase of this product (Rotor-Gene Q, Rotor-Disc) includes a limited, non-transferable license to one or more of US Patents Nos 6,787,338; 7,238,321; 7,081,226; 6,174,670; 6,245,514; 6,569,627; 6,303,305; 6,503,720; 5,871,908; 6,691,041; 7,387,887; and U.S. Patent Applications Nos. 2003-0224434 and 2006-0019253 and all continuations and divisionals, and corresponding claims in patents and patent applications outside the United States, owned by the University of Utah Research Foundation, Idaho Technology, Inc., and/or Roche Diagnostics GmbH, for internal research use or for non-in vitro diagnostics applications. No right is conveyed, expressly, by implication or estoppel, for any reagent or kit, or under any other patent or patent claims owned by the University of Utah Research Foundation, Idaho Technology, Inc., and/or Roche Diagnostics GmbH, or by any other Party. For information on purchasing licences for in-vitro diagnostics applications or reagents, contact Roche Molecular Systems, 4300 Hacienda Drive, Pleasanton, CA 94588, USA.

This real-time thermal cycler (Rotor-Gene Q) is licensed under U.S. Patent No. 6,814,934 and corresponding claims in any foreign counterpart patent thereof owned by Applied Biosystems LLC, in all fields, including research and development, all applied fields, and human and animal in-vitro diagnostics. No rights are conveyed expressly, by implication or estoppel to any patents on real-time methods, including but not limited to 5' nuclease assays, or to any patent claiming a reagent or kit. For further information on purchasing additional rights, contact the Director of Licensing at Applied Biosystems, 850 Lincoln Centre Drive, Foster City, California, 94404, USA.

1056748 01/2010 © 2010 QIAGEN, all rights reserved

www.qiagen.com

Australia = Orders 1-800-243-800 = Fax 03-9840-9888 = Technical 1-800-243-066

Austria = Orders 0800-28-10-10 = Fax 0800-28-10-19 = Technical 0800-28-10-11

Belgium = Orders 0800-79612 = Fax 0800-79611 = Technical 0800-79556

Brazil = Orders 0800-557779 = Fax 55-11-5079-4001 = Technical 0800-557779

Canada = Orders 800-572-9613 = Fax 800-713-5951 = Technical 800-DNA-PREP (800-362-7737)

China = Orders 86-21-3865-3865 = Fax 86-21-3865-3965 = Technical 800-988-0325

Denmark = Orders 80-885945 = Fax 80-885944 = Technical 80-885942

Finland = Orders 0800-914416 = Fax 0800-914415 = Technical 0800-914413

France = Orders 01-60-920-920 = Fax 01-60-920-925 = Technical 01-60-920-930

Germany = Orders 02103-29-12000 = Fax 02103-29-22000 = Technical 02103-29-12400

Hong Kong = Orders 800 933 965 = Fax 800 930 439 = Technical 800 930 425

Ireland = Orders 1800 555 049 = Fax 1800 555 048 = Technical 1800 555 061

Italy = Orders 800-789 544 = Fax 02-334304-826 = Technical 800-787980

Japan = Telephone 03-6890-7300 = Fax 03-5547-0818 = Technical 03-6890-7300

Korea (South) = Orders 080-000-7146 = Fax 02-2626-5703 = Technical 080-000-7145

Luxembourg = Orders 8002-2076 = Fax 8002-2073 = Technical 8002-2067

Mexico = Orders 01-800-7742-639 = Fax 01-800-1122-330 = Technical 01-800-7742-436

The Netherlands = Orders 0800-0229592 = Fax 0800-0229593 = Technical 0800-0229602

Norway = Orders 800-18859 = Fax 800-18817 = Technical 800-18712

Singapore = Orders 1800-742-4362 = Fax 65-6854-8184 = Technical 1800-742-4368

Spain = Orders 91-630-7050 = Fax 91-630-5145 = Technical 91-630-7050

Sweden = Orders 020-790282 = Fax 020-790582 = Technical 020-798328

Switzerland = Orders 055-254-22-11 = Fax 055-254-22-13 = Technical 055-254-22-12

UK = Orders 01293-422-911 = Fax 01293-422-922 = Technical 01293-422-999

USA = Orders 800-426-8157 = Fax 800-718-2056 = Technical 800-DNA-PREP (800-362-7737)

