
PAXgene® Saliva Collector

Do not unscrew the funnel until the saliva sample collection is completed.
Please read the instructions carefully before collecting saliva sample.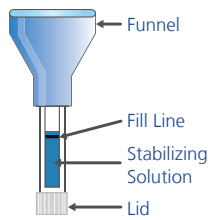

For collection, stabilization, and transportation of 2 ml human saliva.

The PAXgene Saliva Collector is intended for molecular biology applications. This product is not intended for the diagnosis, prevention or treatment of a disease.

Mat. No. 1124399

Materials

The blister package contains one collection device for 2 ml saliva and instructions for use. The device consists of a collection funnel attached to a collection tube and a lid. The collection tube is prefilled with stabilizing solution.

Storage and Shelf-life

Store the PAXgene Saliva Collector at room temperature (15–30°C). When stored properly, the device is stable until the expiration date on the blister packaging label.

Transportation of samples

After the sample is collected, discard the funnel in plastic waste. Close the tube tightly and place it back in the blister packaging. Close the blister flap securely for shipment to return address.

Important Notes

- Make sure that the packaging is intact. If not, do not use the device. Never reuse a device.
- Do not ingest the stabilizing solution. If skin comes in contact with stabilizing solution, wash area with soap and water. If eyes come into contact with stabilizing solution, rinse with water for at least 15 min and consult a physician.
- Rinse mouth 30 min prior collection to remove any food particles in sample.
- Do not drink, eat, smoke, chew gum or brush teeth within 30 min prior to sample collection.
- Assist children and the elderly in using this device.

For an instructional video, please visit www.preanalytix.com

PreAnalytiX GmbH, Feldbachstrasse, CH-8634 Hombrechtikon, Tel.: +49-2103-29-0
Produced in China by Xiamen Zeesan Biotech Co., Ltd. on behalf of PreAnalytiX
(Modified from model SAL-2000L)
Building 1#, No.3701 North Xiang'an Road, (Xiang An) Industrial Area, Torch High-Tech Zone, Xiamen City, Fujian Province, 361101, P.R. China

<p>Do not drink or eat within 30 min prior to sample collection.</p>	<p>1</p> <p>Spit gently into funnel. Saliva collection will take about 5–10 min.</p>
<p>2</p> <p>Bubbles</p> <p>Keep spitting until liquid saliva (excluding bubbles) fills up outer part of the tube and reaches fill line.</p> <p>Shake tube gently to assist saliva flow.</p>	<p>3</p> <p>Hold tube upright and slowly unscrew funnel. This releases stabilizing solution from inner part of the tube. Shake carefully to ensure fluids transfer completely into tube. Discard funnel.</p>
<p>4</p> <p>Keep tube in upright, remove lid from the bottom of tube, place over opening and screw lid on tightly to close.</p>	<p>5</p> <p>Ensure tube is sealed, then invert tube 5 times to mix saliva and stabilizing solution.</p>

For up-to-date licensing information and product-specific disclaimers, see the respective PreAnalytiX® or QIAGEN® kit handbook or user manual. PreAnalytiX and QIAGEN kit handbooks and user manuals are available at www.preanalytix.com and www.qiagen.com or can be requested from PreAnalytiX Technical Services.
HB-2919-001 1124320 08/2021 © 2021 PreAnalytiX GmbH, all rights reserved.

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HB-2919-001 1124320 08/2021 © 2021 PreAnalytiX GmbH, todos los derechos reservados.
HB-2919-001 1124320 08/2021 © 2021 PreAnalytiX GmbH, alle Rechte vorbehalten.