August 2021

# EZ2<sup>®</sup> Connect and EZ2 Connect Fx Safety Instructions Quick-Start Guide

---

# Contents

Introduction.....	3
Safety Information.....	4
Proper use .....	6
Electrical safety .....	12
Environment.....	16
Biological safety.....	19
Chemicals .....	23
Toxic fumes .....	26
Waste disposal .....	27
Mechanical hazards.....	27
Heat hazard .....	29
Radiation safety .....	30
Maintenance safety .....	31
Symbols on the EZ2 .....	36
Installation Procedures.....	42
Site requirements.....	42
Setting up the EZ2 .....	42
To set up the EZ2 .....	43
Connecting the EZ2 via Wi-Fi.....	45
Connecting the EZ2 via LAN cable.....	45
Revision History.....	46

---

# Introduction

The EZ2 Connect and EZ2 Connect Fx (hereafter also referred to as EZ2) is designed to perform automated isolation and purification of nucleic acids in molecular biology applications. The system is intended for users trained in molecular biological techniques and the operation of EZ2.

The EZ2 is intended to be used only in combination with QIAGEN kits indicated for use with the EZ2 Connect or EZ2 Connect Fx.

# Safety Information

Before using the EZ2 , it is essential that you read this guide carefully and pay attention to the safety information. The instructions and safety information in this guide must be followed to ensure safe operation of the instrument and to maintain the instrument in a safe condition.

**Note:** This Quick Start Guide is supplementary to, not a replacement for, the Instrument User Manual

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The advice given in this guide is intended to supplement, not supersede, the normal safety requirements prevailing in the user's country.

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If hazardous material is spilled on or inside the EZ2, the user is responsible for carrying out appropriate decontamination.

**Note:** Do not place items on top of the EZ2.

**Note:** In case of emergency, power OFF the EZ2 and unplug the power cord from the power outlet.

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## Electrical safety

**Note:** If operation of the instrument is interrupted in any way (e.g., due to interruption of the power supply or a mechanical error), first switch off the EZ2 then disconnect the electrical cord from the power supply before attempting troubleshooting or service activity.

<p><b>WARNING</b></p> 	<p><b>Electrical hazard</b></p> <p>Any interruption of the protective conductor (earth/ground lead) inside or outside the instrument or disconnection of the protective conductor terminal is likely to make the instrument dangerous.</p> <p>Intentional interruption is prohibited.</p> <p><b>Lethal voltages inside the instrument</b></p> <p>When the instrument is connected to line power, terminals may be live and opening covers or removing parts is likely to expose live parts.</p>
<p><b>DE</b></p>	<p><b>Stromschlaggefahr</b></p> <p>Jede Unterbrechung des Schutzleiters (Erdungs- bzw. Masseleiter) im Gerät oder außerhalb des Geräts und jede Abtrennung des Schutzleiters am Anschluss der Netzleitung erhöht die Gefahr eines Stromschlags.</p> <p>Eine absichtliche Unterbrechung der Schutzleiterverbindung ist verboten.</p> <p><b>Gefährliche Spannung im Gerät</b></p> <p>Wenn das Gerät an die Stromversorgung angeschlossen ist, sind die Anschlussstellen spannungsführend. Durch das Öffnen der Abdeckungen oder das Entfernen von Gehäuseteilen können spannungsführende Komponenten freigelegt werden.</p>

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To ensure satisfactory and safe operation of the EZ2, follow these guidelines:

- The line power cord must be connected to a line power outlet that has a protective conductor (earth/ground).

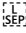
- Place instrument in a location so that the power cord is accessible and can be connected/disconnected.
- Use only the power cord delivered by QIAGEN.
- Do not adjust or replace internal parts of the instrument.
- Do not operate the instrument with any covers or parts removed.
- If liquid has spilled inside the instrument, power OFF the instrument, disconnect it from the power outlet, and contact QIAGEN Technical Services.

If the EZ2 becomes electrically unsafe, prevent other personnel from operating it and contact QIAGEN Technical Services.

The instrument may be electrically unsafe when:

- The EZ2 or the line power cord appears to be damaged.
- The EZ2 has been stored under unfavorable conditions for a prolonged period.
- The EZ2 has been subjected to severe transport stresses.
- Liquids come in contact directly with electrical components of the EZ2.
- The power cord has been exchanged with a non-official power cord.

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<p><b>CAUTION</b></p> 	<p><b>Damage to the instrument</b></p> <p>Direct sunlight may bleach parts of the instrument, cause damage to plastic parts, or interfere with the correct function of the load check. The EZ2 must be located out of direct sunlight.</p>
---	--

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## Biological safety

Specimens and reagents containing materials from humans should be treated as potentially infectious. Use safe laboratory procedures as outlined in publications such as Biosafety in Microbiological and Biomedical Laboratories, HHS (<https://www.cdc.gov/labs/pdf/CDC-BiosafetymicrobiologicalBiomedicalLaboratories-2009-P.pdf>).

Samples may contain infectious agents. You should be aware of the health hazard presented by such agents and should use, store, and dispose of such samples according to the required safety regulations.

**WARNING****Samples containing infectious agents**

Some samples used with the EZ2 may contain infectious agents. Handle such samples with the greatest of care and in accordance with the required safety regulations.

Always wear safety glasses, gloves, and a lab coat.

The responsible persons (e.g., a laboratory manager) must take the necessary precautions to ensure that the surrounding workplace is safe and that the instrument operators are suitably trained and not exposed to hazardous levels of infectious agents as defined in the applicable Safety Data Sheets (SDSs) or the OSHA\*, ACGIH†, or COSHH‡ documents.

Venting for fumes and disposal of wastes must be in accordance with all national, state, and local health and safety regulations and laws.

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\* OSHA: Occupational Safety and Health Administration (United States of America).

† ACGIH: American Conference of Government Industrial Hygienists (United States of America).

‡ COSHH: Control of Substances Hazardous to Health (United Kingdom).

## Chemicals

<p><b>WARNING</b></p> 	<p><b>Hazardous chemicals</b></p> <p>Some chemicals used with the EZ2 may be hazardous or may become hazardous after completion of a purification.</p> <p>Always wear safety glasses, gloves, and a lab coat.</p> <p>The responsible persons (e.g., a laboratory manager) must take the necessary precautions to ensure that the surrounding workplace is safe and that the instrument operators are not exposed to hazardous levels of toxic substances (chemical or biological) as defined in the applicable Safety Data Sheets (SDSs) or the OSHA*, ACGIH†, or COSHH‡ documents.</p> <p>Venting for fumes and disposal of wastes must be in accordance with all national, state, and local health and safety regulations and laws.</p>
---	---

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**Note:** If you work with volatile solvents, toxic substances, etc., you must provide an efficient laboratory ventilation system to remove vapors that may be produced.

## Waste disposal

Used consumables, such as reagent cartridges and disposable filter-tips, may contain hazardous chemicals or infectious agents from the purification process. Such waste must be collected and disposed of properly, according to local safety regulations.

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## Mechanical hazards

The hood of the EZ2 must remain closed during operation of the instrument. Only open the hood when instructed to do so by the instructions for use.

The worktable of the EZ2 moves during operation of the instrument. While loading the worktable, always stand clear of the instrument. Do not lean on the worktable when the robotic arm of the instrument is moving to reach loading position with its lid open. Wait until the robotic arm completed its movements before you start to load or unload.

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## Heat hazard

The EZ2 worktable contains a heating system.

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## Radiation safety

The EZ2 has a UV LED lamp. The wavelength of the UV light produced by the UV LED lamp is 270 to 285 nm. This wavelength corresponds to ultraviolet light type C, which can be used for decontamination procedures. A mechanical lock ensures that the hood must be closed for operation of the UV LED. If the hood sensor or lock is not functioning correctly, contact QIAGEN Technical Services.

<b>WARNING</b> 	<b>UV radiation</b> Avoid looking directly into UV light. Do not expose your skin to UV light.
<b>DE</b>	<b>UV-Strahlung</b> Vermeiden Sie es, direkt in das UV-Licht zu schauen. Setzen Sie Ihre Haut nicht dem UV-Licht aus.
<b>FR</b>	<b>Rayonnement UV</b> Éviter de regarder directement la lumière UV. Ne pas exposer votre peau à la lumière UV.

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# Symbols on the EZ2

The following symbols appear on the EZ2.

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---

# Installation Procedures

This section provides you with important information on how to install the EZ2, including the required instrument setup and network settings. For detailed unpacking and installing information, refer to the respective user manual.

The EZ2 is a plug-and-play instrument. Unpacking and installation procedures are easy to follow, but a person who is familiar with laboratory equipment should oversee the installation.

## Site requirements

The EZ2 must be located out of direct sunlight, away from heat sources, and away from sources of vibration and electrical interference. Refer to the *EZ2 Connect and EZ2 Connect Fx User Manual* for the operating conditions (temperature and humidity) and detailed site, power, and grounding requirements.

## Setting up the EZ2

This section describes important actions that must be performed before operating the EZ2.

The following items are delivered with the EZ2 Connect and EZ2 Connect Fx:

- EZ2 Connect or EZ2 Connect Fx instrument
- Quick-start guide(this document)
- Left and right cartridge racks
- Left and right tip racks
- Power cord set
- USB drive
- Silicone grease

---

Additional items are delivered with the EZ2 Connect Fx:

- Handheld barcode scanner
- Left and right tip racks for flip cap tubes

To set up the EZ2

1. Before unpacking the EZ2, move the package to the site of installation and check that the arrows on the package point upward. In addition, check whether the package is damaged. In case of damage, contact QIAGEN Technical Services.
2. Open the top of the transportation box to remove the top layer (PE foam).
3. Remove the accessories box together with the PE foam at its perimeter.
4. Remove the outer carton box by holding at the cut out area and lift up the outer carton box.
5. Remove two pieces of protector from the unit.
6. Move the unit to the workbench or trolley from the packaging. When lifting the EZ2, slide your fingers under the side of the instrument and keep your back straight.

**Important:** Two persons are required to lift the EZ2.

**Important:** Do not hold the touchscreen display while unpacking or lifting the EZ2, this might damage the instrument.

7. Remove the transport lock for P-axis by pulling the transport lock from bottom.
8. Remove silica gel from the unit.
9. Remove the transport lock for Y-axis by pushing the bottom of transport lock toward rear direction and pull it out from the rear. There is a total of two transport lock for Y-axis.
10. Check if the packing list document is included after unpacking the EZ2.
11. Read the packing list to check that you have received all items. If anything is missing, contact QIAGEN Technical Services.

- 
12. Check that the EZ2 is not damaged and that there are no loose parts. If anything is damaged, contact QIAGEN Technical Services. Make sure that the EZ2 has equilibrated to ambient temperature before operating it.
  13. Retain the package in case you need to transport the EZ2 in the future. Refer to section “Packing the EZ2” of the user manual for more details. Using the original package minimizes the possibility of damage during transportation of the EZ2.
  14. Before turning on the instrument, plug the Wi-Fi adapter into one of the USB ports located on the front of the instrument or on the rear of the touchscreen.
  15. Connect the barcode scanner to one of the USB ports located on the front of the instrument or on the rear of the touchscreen.
  16. Connect the EZ2 to a power outlet using the provided power cord.
  17. To power on the EZ2, first ensure that the hood is closed then press the power button. The startup screen appears, and the instrument is initialized. After the initialization is done, the Login screen appears.
  18. To log in for the first time, enter **Admin** in both **User ID** and **Password** fields. After this login, you will have administrator rights with the option to set up other users. Refer to the instrument user manual for more details on user management.
  19. From the **Configuration** menu under the **System** tab, you can modify the **Device Name**, **Date**, and **Time** fields.
  20. You may adjust the settings of the EZ2 according to your preferences through the **Configuration** menu, **Device Settings** tab.
  21. Before first use in the final installation site, the camera exposure calibration **must** be performed through the **Maintenance** menu, **Camera LED** tab (EZ2 Connect Fx only).

---

## Connecting the EZ2 via Wi-Fi

The EZ2 can be connected to a Wi-Fi network using the Wi-Fi adapter inserted into one of the USB ports. The adapter supports the 802.11b, 802.11g, and 802.11n Wi-Fi standards, and WEP, WPA-PSK, and WPA2-PSK encryption.

- From the **Configuration** menu under the **Wi-Fi** tab, you can set up the Wi-Fi network.
- Scan for available networks.
- From the list, select one of the available networks.
- Enter the Wi-Fi password and connect.

## Connecting the EZ2 via LAN cable

From the **Configuration** menu under the **LAN** tab, enter the local network settings to set up the LAN network.

**Note:** You might need the help of an IT professional to enter detailed network settings.

# Revision History

Revision	Description
08/2021	First version of the Quick-Start Guide.

For further support with technical difficulties, please contact QIAGEN Technical Services at [www.qiagen.com](http://www.qiagen.com).

For up-to-date licensing information and product-specific disclaimers, see the respective QIAGEN user manual. QIAGEN kit handbooks and user manuals are available at [www.qiagen.com](http://www.qiagen.com) or can be requested from QIAGEN Technical Services or your local distributor.

---

## Notes

---

Trademarks: QIAGEN®, Sample to Insight®, EZ2® (QIAGEN Group). Registered names, trademarks, etc. used in this document, even when not specifically marked as such, are not to be considered unprotected by law.

HB-2925-001 08/2021 1125025 © 2021 QIAGEN, all rights reserved.