

November 2018

miScript[®] miRNA PCR Array Handbook

miScript II RT Kit

miScript SYBR[®] Green PCR Kit

miScript miRNA PCR Arrays

miScript miRNA QC PCR Array

For SYBR Green-based, real-time PCR

profiling of microRNAs using pathway-focused
arrays, HC arrays, and miRNome arrays

Contents

Kit Contents.....	4
Storage	8
Intended Use	9
Safety Information.....	9
Quality Control.....	9
Introduction	10
Principle and Procedure.....	14
Template RNA requirements.....	26
Equipment and Reagents to Be Supplied by User	27
Protocol: Reverse Transcription for Quantitative, Real-Time PCR	29
Protocol: Real-Time PCR for Mature miRNA Expression Profiling	33
Protocol: Data Analysis.....	40
Protocol: cDNA Quality Control Prior to Profiling Mature miRNA.....	43
Protocol: Data Analysis for Quality Control Using miScript miRNA QC PCR Arrays	46
Troubleshooting Guide	49
Appendix A: Real-Time PCR Data Output and Dissociation Curve Analysis	52
Appendix B: General Remarks on Handling RNA	56
Handling RNA	56
General handling	56
Disposable plasticware	57
Glassware	57
Solutions	57

Appendix C: Preparation, Quantification, and Storage of RNA.....	59
RNA preparation and quality.....	59
Determining concentration and purity of RNA.....	59
Storage of RNA.....	60
Ordering Information	61

Kit Contents

miScript II RT Kit	(12)	(50)
Catalog no.	218160	218161
Number of standard reactions*	12	50
miScript Reverse Transcriptase Mix	24 µl	100 µl
10x miScript Nucleics Mix	50 µl	200 µl
5x miScript HiSpec Buffer	100 µl	400 µl
5x miScript HiFlex Buffer	100 µl	400 µl
RNase-Free Water	1.9 ml	1.9 ml
Quick-Start Protocol	1	1

* A standard reaction is 20 µl in volume with up to 2 µg total RNA (when using miScript HiSpec Buffer).

miScript SYBR Green PCR Kit	(200)	(1000)
Catalog no.	218073	218075
Number of 50 µl reactions	200[†]	1000[‡]
2x QuantiTect® SYBR Green PCR Master Mix, containing: HotStarTaq® DNA Polymerase QuantiTect SYBR Green PCR Buffer dNTP mix, including dUTP SYBR Green I ROX™ passive reference dye 5 mM MgCl ₂	3 x 1.7 ml	25 ml
10x miScript Universal Primer	1 ml	5 x 1 ml
RNase-Free Water	2 x 2 ml	20 ml
Quick-Start Protocol	2	2

[†] Provides sufficient reagents for 4 x 96-well arrays (110 µl residual), 2 x 384-well pathway-focused arrays (1210 µl residual), 2 x 384-well HC/miRNome arrays (1510 µl residual), or 5 Rotor-Disc 100 arrays (110 µl residual).

[‡] Provides sufficient reagents for 20 x 96-well arrays (0 µl residual), 12 x 384-well pathway-focused arrays (1100 µl residual), 13 x 384-well HC/miRNome arrays (850 µl residual), or 25 Rotor-Disc 100 arrays (0 µl residual).

Pathway-Focused miScript miRNA PCR Array							
Catalog no. Format	Varies A	C	D	E	F	G	R
96-well plate containing dried assays	2, 12, or 24	2, 12, or 24	2, 12, or 24	–	2, 12, or 24	–	–
384-well plate containing dried assays	–	–	–	4	–	4	–
Rotor-Disc® 100 containing dried assays	–	–	–	–	–	–	2, 12, or 24
Optical Thin-Wall 8-Cap Strips (12 per plate)	24, 144, or 288	–	24, 144, or 288	–	–	–	–
Optical Adhesive Film (1 per plate)	–	2, 12, or 24	–	4	2, 12, or 24	4	–
384EZLoad Covers (1 set of 4 per plate)	–	–	–	4 sets	–	4 sets	–
Rotor-Disc Heat Sealing Film (1 per Rotor-Disc)	–	–	–	–	–	–	2, 12, or 24

miScript miRNA HC PCR Array			
Catalog no. Format	Varies	E	G
384-well plate containing dried assays		2, 12, or 24	2, 12, or 24
Optical Adhesive Film (1 per plate)		2, 12, or 24	2, 12, or 24

miRNome miScript miRNA PCR Array							
Catalog no. Format	Varies A	C	D	E	F	G	R
96-well plate containing dried assays	Varies	Varies	Varies	–	Varies	–	–
384-well plate containing dried assays	–	–	–	Varies	–	Varies	–
Rotor-Disc 100 containing dried assays	–	–	–	–	–	–	Varies
Optical Thin-Wall 8-Cap Strips (12 per plate)	Varies	–	Varies	–	–	–	–
Optical Adhesive Film (1 per plate)	–	Varies	–	Varies	Varies	Varies	–
Rotor-Disc Heat Sealing Film (1 per Rotor-Disc)	–	–	–	–	–	–	Varies

miScript miRNA QC PCR Array							
Catalog no. Format	Varies A	C	D	E	F	G	R
96-well plate containing dried assays	1	1	1	-	1	-	-
384-well plate containing dried assays	-	-	-	1	-	1	-
Rotor-Disc 100 containing dried assays	-	-	-	-	-	-	1
Optical Thin- Wall 8-Cap Strips	12	-	12	-	-	-	-
Optical Adhesive Film	-	1	-	1	1	1	-
Rotor-Disc Heat Sealing Film (1 per Rotor-Disc)	-	-	-	-	-	-	1

Cyclers for use with array formats

Format	Suitable real-time cyclers	Plate
A	Applied Biosystems® models 5700, 7000, 7300, 7500, 7700, 7900HT, ViiA™ 7 (96-well block); Bio-Rad® models iCycler®, iQ™5, MyiQ™, MyiQ2; Bio-Rad/MJ Research Chromo4™; Eppendorf® Mastercycler® ep realplex models 2, 2S, 4, 4S; Stratagene® models Mx3005P®, Mx3000P®; Takara TP-800	96-well
C	Applied Biosystems models 7500 (Fast block), 7900HT (Fast block), StepOnePlus™, ViiA 7 (Fast block)	96-well
D	Bio-Rad CFX96™; Bio-Rad/MJ Research models DNA Engine Opticon®, DNA Engine Opticon 2; Stratagene Mx4000®	96-well
E	Applied Biosystems models 7900HT (384-well block), ViiA 7 (384-well block); Bio-Rad CFX384™	384-well
F	Roche® LightCycler® 480 (96-well block)	96-well
G	Roche LightCycler 480 (384-well block)	384-well
R	Rotor-Gene® Q; Rotor-Gene 6000; other Rotor-Gene cyclers	Rotor-Disc 100

Storage

The miScript II RT Kit and miScript SYBR Green PCR Kit are shipped on dry ice. The kits, including all reagents and buffers, should be stored immediately upon receipt at -20°C in a constant-temperature freezer.

miScript miRNA PCR Arrays and miScript miRNA QC PCR Arrays are shipped at ambient temperature, on ice, or on dry ice depending on the destination and accompanying products. Upon receipt, store at -20°C . If stored under these conditions, miScript miRNA PCR Arrays and miScript miRNA QC PCR Arrays are stable for 6 months after receipt.

Intended Use

The miScript II RT Kit, miScript SYBR Green PCR Kit, miScript miRNA PCR Arrays and miScript miRNA QC PCR Array are intended for molecular biology applications. These products are not intended for the diagnosis, prevention or treatment of a disease.

All due care and attention should be exercised in the handling of the products. We recommend all users of QIAGEN® products to adhere to the NIH guidelines that have been developed for recombinant DNA experiments, or to other applicable guidelines.

Safety Information

When working with chemicals, always wear a suitable lab coat, disposable gloves and protective goggles. For more information, please consult the appropriate safety data sheets (SDSs). These are available online in convenient and compact PDF format at www.qiagen.com/us/support/qa-qc-safety-data/safety-data-sheets/ where you can find, view and print the SDS for each QIAGEN kit and kit component.

Quality Control

In accordance with QIAGEN's ISO-certified Quality Management System, each lot of miScript II RT Kit and miScript SYBR Green PCR Kit is tested against predetermined specifications to ensure consistent product quality.

Introduction

The miScript PCR System consists of the miScript II RT Kit, miScript PreAMP PCR Kit, miScript miRNA PCR Array, miScript Primer Assay, miScript SYBR Green PCR Kit, and miScript miRNA PCR Array data analysis tool. The miScript PCR System allows sensitive and specific detection and quantification of microRNA (miRNA). The miScript PCR System uses total RNA that contains miRNA as the starting material for cDNA synthesis, and separate enrichment of small RNA is not needed. A single cDNA preparation can be used with a miScript miRNA PCR Array to rapidly profile the expression of mature miRNAs.

miScript miRNA PCR Arrays are mature miRNA-specific forward primers (miScript Primer Assays) that have been arrayed in biologically relevant pathway-focused and whole miRNome panels. These PCR arrays are provided in ready-to-use 384-well plate, 96-well plate, and 100-well Rotor-Disc formats. miScript miRNA PCR Arrays, which are available for several species, provide guaranteed high performance and are fully customizable. Each assay in a miScript miRNA PCR Array has been verified to ensure sensitive and specific detection of mature miRNA by real-time PCR. A free, Web-based miScript miRNA PCR Array data analysis tool simplifies the analysis of real-time PCR data. Once raw threshold cycle (C_T) data has been uploaded, the tool automatically performs all fold-change calculations using the $\Delta\Delta C_T$ method of relative quantification and presents the results in several formats. Mature miRNome expression profiling is now within reach of every laboratory because of the ease, convenience, and consistent performance of miScript miRNA PCR Arrays. miScript miRNA PCR Arrays are at the forefront of real-time PCR-based mature miRNA profiling tools.

miScript miRNA PCR Array workflow

Prepare reverse-transcription reaction




Incubate at 37°C for 60 min,
then at 95°C for 5 min


Prepare PCR mix






Add PCR mix to miScript miRNA PCR Array

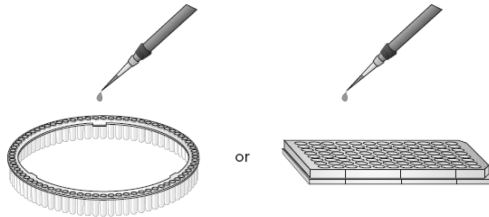

Perform real-time PCR

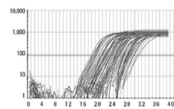

Analyze results using miScript miRNA PCR Array data analysis tool

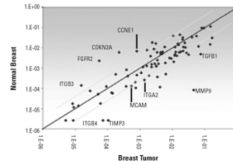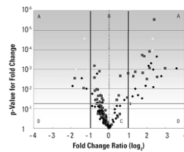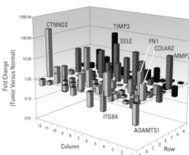

miScript II RT Kit

The expanded miScript II RT Kit includes miScript Reverse Transcriptase Mix, 10x miScript Nucleics Mix, 5x miScript HiSpec Buffer, and 5x miScript HiFlex Buffer. miScript Reverse Transcriptase Mix is an optimized blend of poly(A) polymerase and reverse transcriptase. 10x miScript Nucleics Mix contains dNTPs, rATP, oligo-dT primers, and an internal synthetic RNA control (miRNA reverse transcription control [miRTC]) that is used to assess reverse transcription performance during profiling experiments with miScript miRNA PCR Arrays. Two buffers, 5x miScript HiSpec Buffer and 5x miScript HiFlex Buffer, are provided in the miScript II RT Kit to meet the distinctive needs of miRNA quantification studies using real-time PCR (Figure 1). The unique, patent-pending formulation of 5x miScript HiSpec Buffer facilitates the selective conversion of mature miRNAs into cDNA, which can then be used for miRNA quantification with either miScript miRNA PCR Arrays or miScript Primer Assays. Using miScript HiSpec Buffer, the conversion of long RNAs, such as mRNAs, is suppressed. As a result, background signals potentially contributed by long RNAs are nonexistent, making a reverse-transcription reaction buffered using miScript HiSpec Buffer highly suitable for profiling the expression of mature miRNAs. Only 5x miScript HiSpec Buffer should be used to prepare cDNA for real-time PCR with miScript miRNA PCR Arrays. For protocols using miScript Primer Assays, refer to the *miScript PCR System Handbook*. If working with samples that contain low amounts of RNA, 5x miScript HiSpec Buffer should also be used to prepare cDNA for preamplification with the miScript PreAMP PCR Kit. For protocols using the miScript PreAMP PCR Kit, refer to the *miScript PreAMP Handbook*.

5x miScript HiFlex Buffer promotes conversion of all RNA species (mature miRNA, precursor miRNA, other noncoding RNA, and mRNA) into cDNA, and this cDNA can then be used to quantify each RNA species (using appropriate primer assays). Do not use 5x miScript HiFlex Buffer to prepare cDNA for real-time PCR with miScript miRNA PCR Arrays. For protocols using miScript HiFlex Buffer, refer to the *miScript PCR System Handbook*.

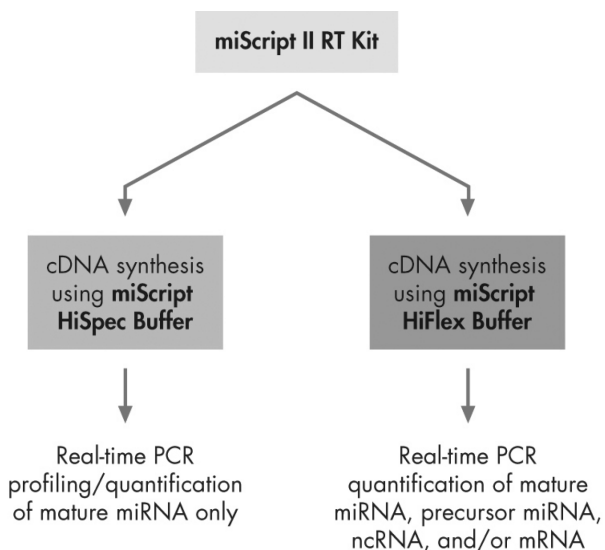


Figure 1. Mature miRNA, precursor miRNA, other noncoding RNA, and mRNA detection.

Two buffers are provided with the miScript II RT Kit. Use miScript HiSpec Buffer for cDNA synthesis to enable either mature miRNA profiling (using miScript miRNA PCR Arrays) or mature miRNA quantification (using individual miScript Primer Assays). For protocols using miScript HiSpec Buffer in combination with miScript Primer Assays, refer to the miScript PCR System Handbook. If samples being assessed have low RNA content, miScript HiSpec Buffer should be used to prepare cDNA for preamplification with the miScript PreAMP PCR Kit. For protocols using the miScript PreAMP PCR Kit, refer to the *miScript PreAMP Handbook*. Use miScript HiFlex Buffer for cDNA synthesis to enable quantification of mature miRNA, precursor miRNA, other noncoding RNA (ncRNA), and/or mRNA from the same cDNA (using appropriate primer assays). For protocols using miScript HiFlex Buffer, refer to the *miScript PCR System Handbook*. Only 5x miScript HiSpec Buffer should be used to prepare cDNA for real-time PCR with miScript miRNA PCR Arrays.

Principle and Procedure

Mature miRNAs are naturally occurring, 22-nucleotide, noncoding RNAs that mediate posttranscriptional gene regulation. Unlike mRNAs, miRNAs are not polyadenylated in nature.

Reverse transcription using miScript HiSpec Buffer

When reverse-transcription reactions are performed in miScript HiSpec Buffer, mature miRNAs and certain small nucleolar RNAs and small nuclear RNAs (snoRNAs and snRNAs, see “Controls in miScript miRNA PCR Arrays and miScript miRNA QC PCR Arrays”) are selectively converted into cDNA. The conversion of long RNAs, such as mRNAs, is suppressed. Mature miRNAs are polyadenylated by poly(A) polymerase and reverse transcribed into cDNA using oligo-dT primers (Figure 2). Polyadenylation and reverse transcription are performed in parallel in the same tube. The oligo-dT primers have a 3' degenerate anchor and a universal tag sequence on the 5' end, allowing amplification of mature miRNA in the real-time PCR step. miScript miRNA PCR Arrays, used in combination with the miScript SYBR Green PCR Kit, enable quantification of mature miRNA by real-time PCR. The combination of polyadenylation and the universal tag addition ensures that miScript miRNA PCR Arrays do not detect genomic DNA.

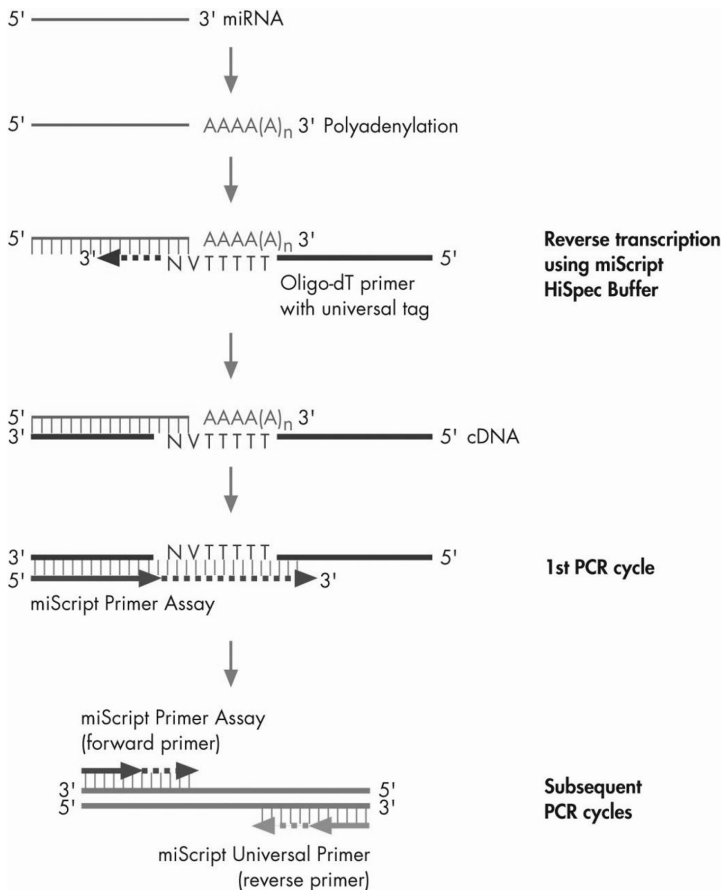


Figure 2. Conversion of mature miRNAs into cDNA and subsequent detection.

In a reverse-transcription reaction using miScript HiSpec Buffer, mature miRNAs are polyadenylated by poly(A) polymerase and subsequently converted into cDNA by reverse transcriptase with oligo-dT priming. The cDNA is then used for real-time PCR profiling of mature miRNA expression (using a miScript miRNA PCR Array and the miScript Universal Primer).

Optional cDNA preamplification using miScript PreAMP PCR Kit

The miScript PreAMP PCR Kit enables researchers to perform miRNA profiling experiments using very limited amounts of starting RNA. This is particularly important when working with samples such as body fluids, formalin-fixed, paraffin-embedded (FFPE) samples, and small cell number samples such as laser capture microdissection (LCM) samples, flow-sorted cells, circulating tumor cells, and fine needle biopsies, where the low RNA yields obtained are often insufficient for reliable miRNA profiling experiments.


The miScript PreAMP PCR Kit, in combination with miScript PreAMP Primer Mixes, uses highly multiplex, PCR-based preamplification of up to 400 miRNA-specific cDNA targets in one reaction. This breakthrough technology enables accurate and comprehensive miRNome-wide expression analysis with as little as 10 ng total RNA. For more information, refer to the *miScript PreAMP Handbook* or visit www.qiagen.com/miRNA.

Mature miRNA expression profiling using miScript miRNA PCR Arrays

cDNA prepared in a reverse-transcription reaction using miScript HiSpec Buffer serves as the template for real-time PCR analysis using a miScript miRNA PCR Array, which contains miRNA-specific miScript Primer Assays, and the miScript SYBR Green Kit, which contains the miScript Universal Primer (reverse primer) and QuantiTect SYBR Green PCR Master Mix. To profile mature miRNA expression, a premix of cDNA, miScript Universal Primer, QuantiTect SYBR Green PCR Master Mix, and RNase-free water is added to a miScript miRNA PCR Array.

miScript miRNA PCR Array plate layout

miScript miRNA PCR Arrays are available in 96-well, 384-well, and Rotor-Disc 100 formats (Figures 3–6). Each array contains several control assays. The purpose of each control is described on page 24.

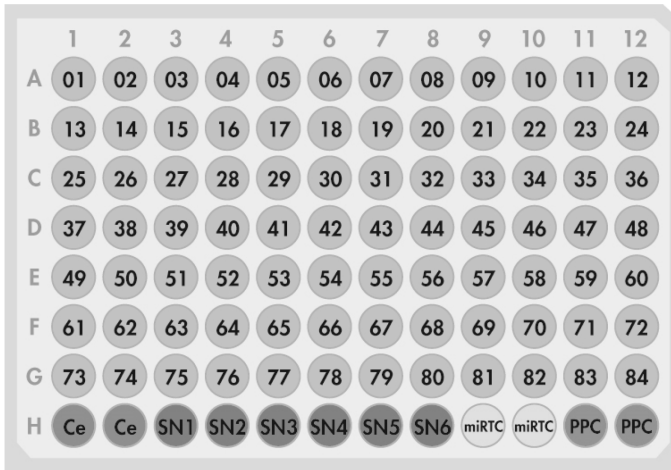

C. elegans miR-39 miScript Primer Assay snoRNA/snRNA miScript PCR Controls Reverse transcription control Positive PCR control

Figure 3. Pathway-Focused or miRNome miScript miRNA PCR Array layout for plate formats A, C, D, F.

Wells A1 to G12 (1–84) each contain a miScript Primer Assay for a pathway/disease/functionally-related mature miRNA. Wells H1 and H2 contain replicate *C. elegans* miR-39 miScript Primer Assays that can be used as an alternative normalizer for array data (Ce). Wells H3 to H8 each contain an assay for a different snoRNA/snRNA that can be used as a normalization control for the array data (SN1=SNORD61 assay, SN2=SNORD68 assay, SN3=SNORD72 assay, SN4=SNORD95 assay, SN5=SNORD96A assay, SN6=RNU6B/RNU6-2 assay). Wells H9 and H10 contain replicate miRTC miScript Primer Assays (miRTC). Wells H11 and H12 contain replicate positive PCR controls (PPC).

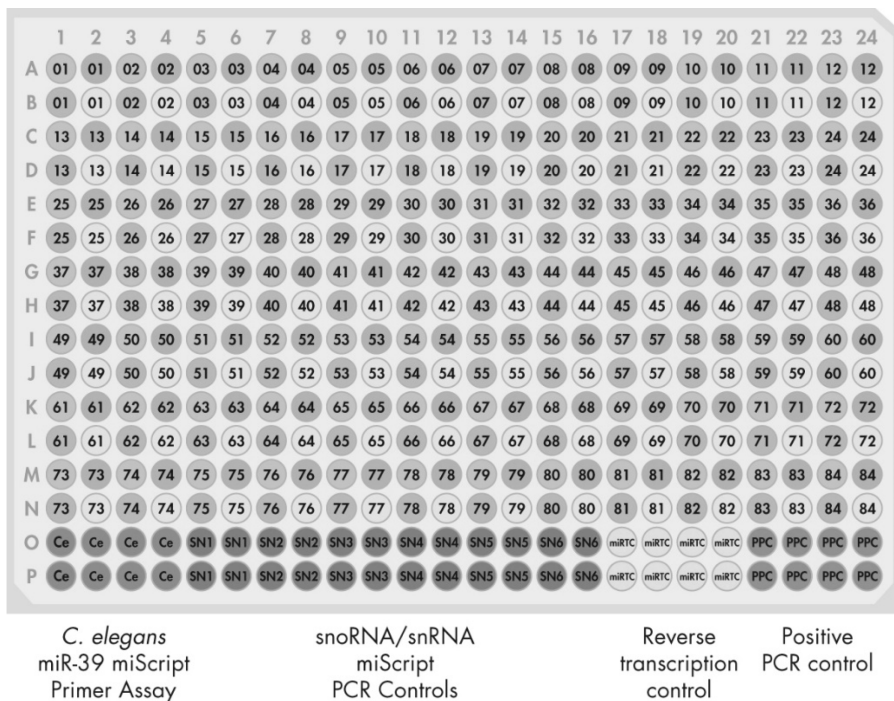


Figure 4. Pathway-Focused miScript miRNA PCR Array layout for plate formats E, G.

Pathway-Focused miScript miRNA PCR Arrays in formats E and G include 4 replicates of the same assays as provided in the 96-well format shown in Figure 3.

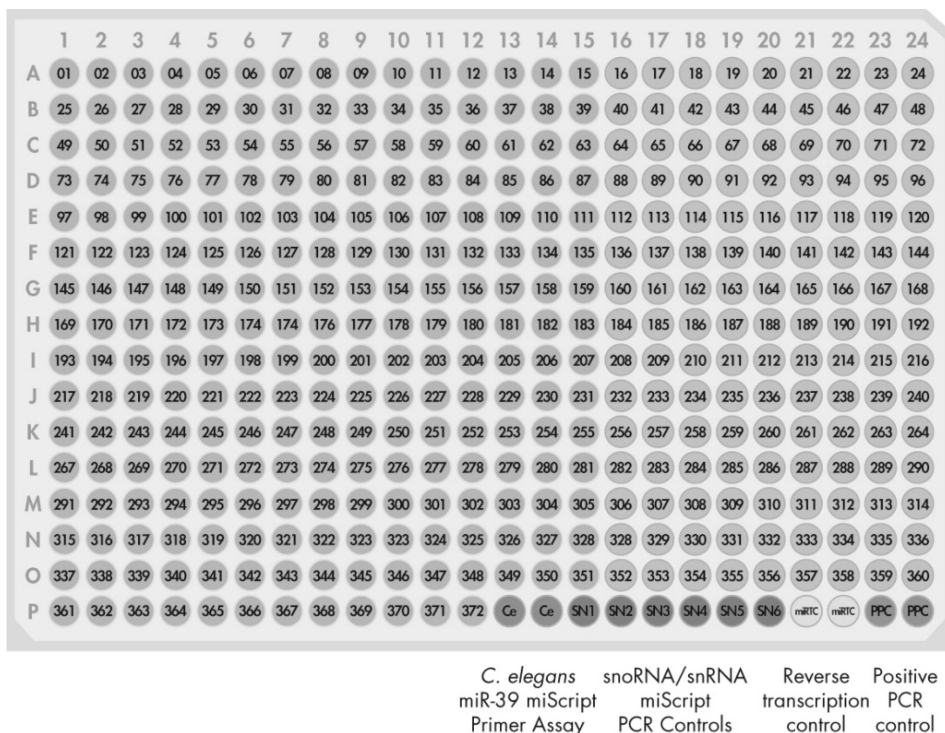


Figure 5. miRNome miScript miRNA PCR Array and miScript miRNA HC PCR Array layout for plate formats E, G. Wells A1 to P12 (1–372) each contain a miScript Primer Assay for a miRNome or pathway/disease/functionally-related mature miRNA. Wells P13 and P14 contain replicate *C. elegans* miR-39 miScript Primer Assays that can be used as an alternative normalizer for array data (Ce). Wells P15 to P20 each contain an assay for a different snoRNA/snRNA that can be used as a normalization control for the array data (SN1=SNORD61 assay, SN2=SNORD68 assay, SN3=SNORD72 assay, SN4=SNORD95 assay, SN5=SNORD96A assay, SN6=RNU6B/RNU6-2 assay). Wells P21 and P22 contain replicate miRTC miScript Primer Assays (miRTC). Wells P23 and P24 contain replicate positive PCR controls (PPC).


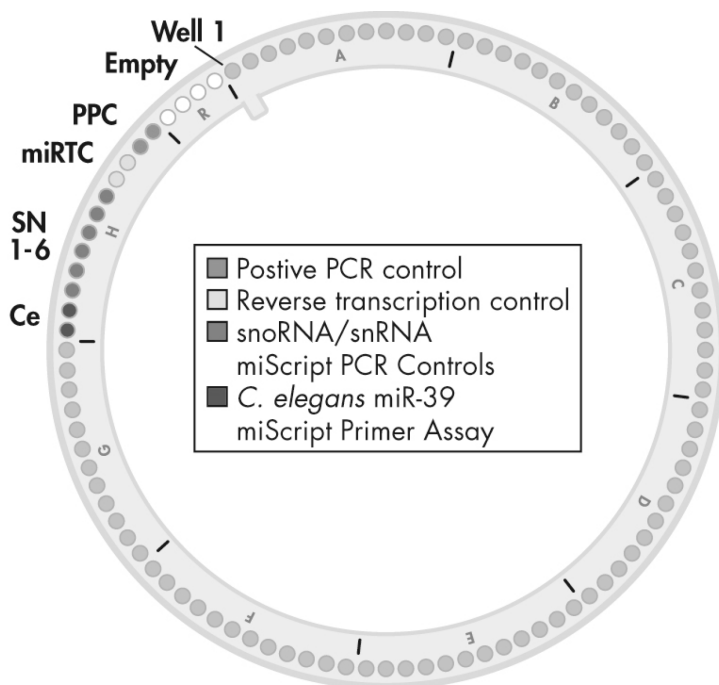


Figure 6. miScript miRNA PCR Array layout for Rotor-Disc format R.

Wells 1 to 84 each contain a miScript Primer Assay for a pathway/disease/functionally-related gene. Wells 85 and 86 contain replicate *C. elegans* miR-39 miScript Primer Assays that can be used as an alternative normalizer for array data (**Ce**). Wells 87 to 92 each contain an assay for a different snoRNA/snRNA that can be used as a normalization control for the array data (**SN1**=SNORD61 assay, **SN2**=SNORD68 assay, **SN3**=SNORD72 assay, **SN4**=SNORD95 assay, **SN5**=SNORD96A assay, **SN6**=RNU6B/RNU6-2 assay). Wells 93 and 94 contain replicate miRTC miScript Primer Assays (**miRTC**). Wells 95 and 96 contain replicate positive PCR controls (**PPC**). Wells 97–100 are empty.

miScript miRNA QC PCR Array plate layout

miScript miRNA QC PCR Arrays enable assessment of the quality of multiple cDNA samples using real-time PCR. miScript miRNA QC PCR Arrays are available in 96-well, 384-well, and Rotor-Disc 100 formats (Figures 7–9). The purpose of each control is described on page 24. The 96-well plate and Rotor-Disc 100 formats allow quality control of up to 8 cDNA samples. The 384-well format allows quality control of up to 32 cDNA samples.

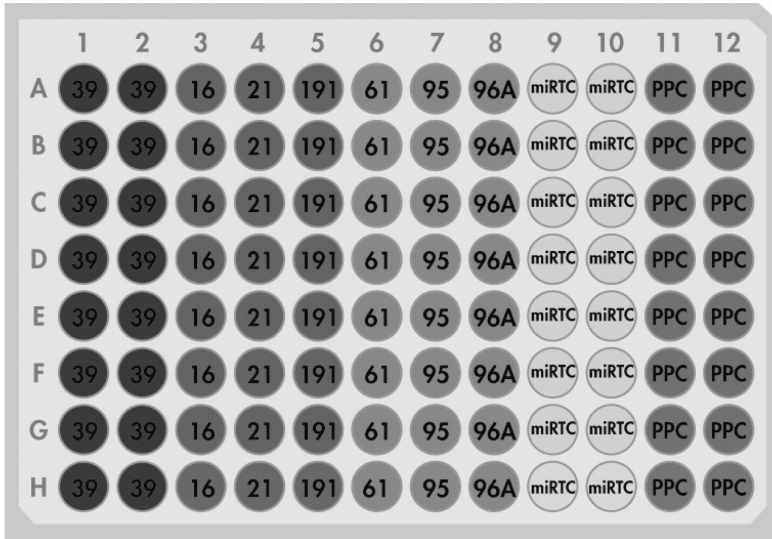

C. elegans miScript Primer Assay snoRNA miScript PCR Controls miRTC miScript Primer Assay Positive PCR control

Figure 7. miScript miRNA QC PCR Array layout for plate formats A, C, D, F.

Wells 1 and 2 of each row contain replicate *C. elegans* miR-39 miScript Primer Assays (39). Well 3 of each row contains a miR-16 miScript Primer Assay (16). Well 4 of each row contains a miR-21 miScript Primer Assay (21). Well 5 of each row contains a miR-191 miScript Primer Assay (191). Wells 6 to 8 of each row each contain an assay for a different snoRNA (61 = SNORD61 assay, 95 = SNORD95 assay, 96A = SNORD96A assay). Wells 9 and 10 of each row contain replicate miRTC miScript Primer Assays (miRTC). Wells 11 and 12 of each row contain replicate positive PCR controls (PPC). These formats enable the quality assessment of up to 8 cDNA samples.

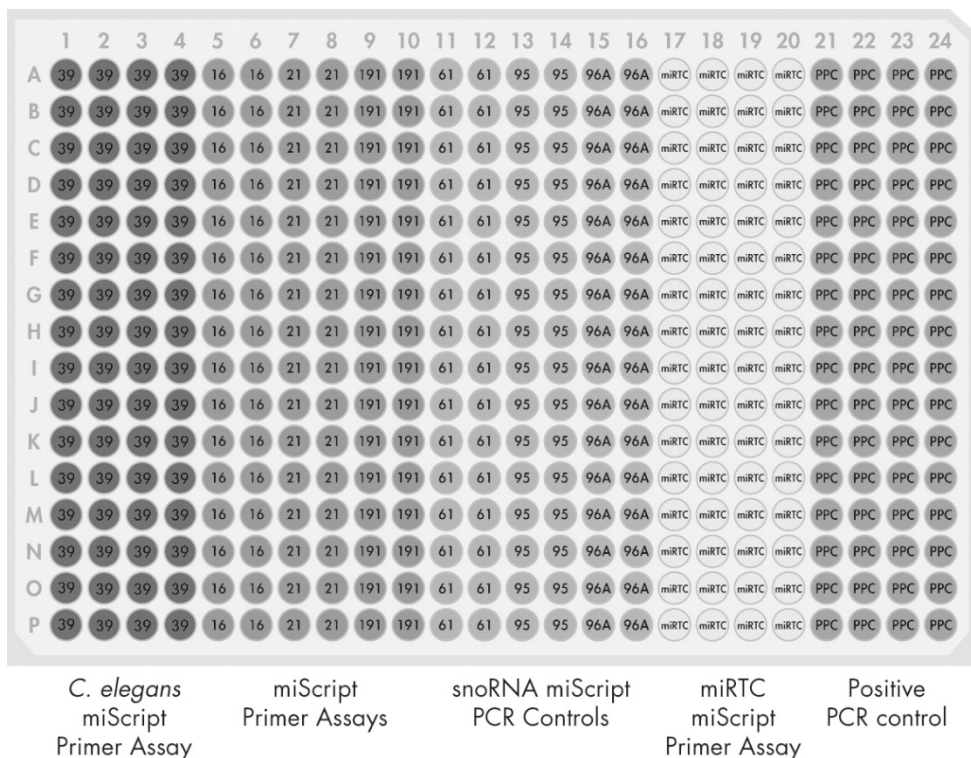


Figure 8. miScript miRNA QC PCR Array layout for plate formats E and G.

Wells 1 to 4 of each row contain replicate *C. elegans* miR-39 miScript Primer Assays (39). Wells 5 and 6 of each row contain replicate miR-16 miScript Primer Assays (16). Wells 7 and 8 of each row contain replicate miR-21 miScript Primer Assays (21). Wells 9 and 10 of each row contain replicate miR-191 miScript Primer Assays (191). Wells 11 to 16 of each row contain assays for different snoRNAs (61 = SNORD61 assay, 95 = SNORD95 assay, 96A = SNORD96A assay). Wells 17 to 20 of each row contain replicate miRTC miScript Primer Assays (miRTC). Wells 21 to 24 of each row contain replicate positive PCR controls (PPC). These formats enable the quality assessment of up to 32 cDNA samples.

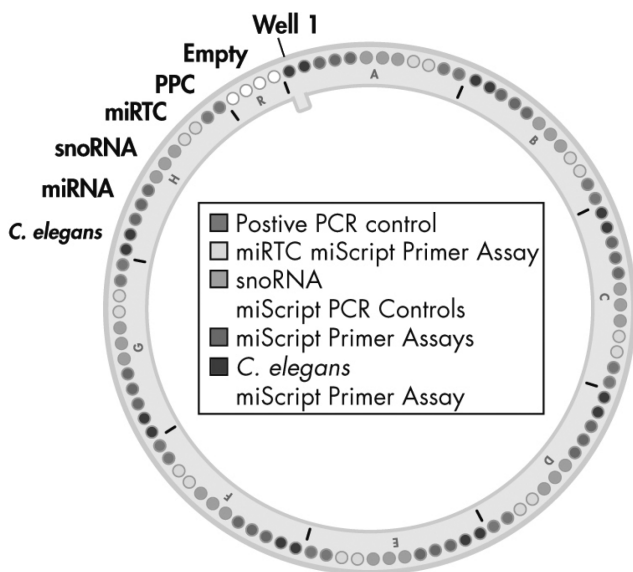


Figure 9. miScript miRNA QC PCR Array layout for Rotor-Disc format R.

Wells 1 and 2 contain replicate *C. elegans* miR-39 miScript Primer Assays. Well 3 contains a miR-16 miScript Primer Assay. Well 4 contains a miR-21 miScript Primer Assay. Well 5 contains a miR-191 miScript Primer Assay. Wells 6 to 8 each contain an assay for a different snoRNA (Well 6 = SNORD61 assay, Well 7 = SNORD95 assay, Well 8 = SNORD96A assay). Wells 9 and 10 contain replicate miRTC miScript Primer Assays (miRTC). Wells 11 and 12 contain replicate positive PCR controls (PPC). This pattern is repeated 7 additional times from wells 13 to 96. Wells 97 to 100 are empty. This format enables the quality assessment of up to 8 cDNA samples.

Controls in miScript miRNA PCR Arrays and miScript miRNA QC PCR Arrays

The final 12 wells of each miScript miRNA PCR Array contain controls. The purpose of the controls in miScript miRNA PCR Arrays and the miScript miRNA QC PCR Array is detailed below and in Table 1.

Table 1. Controls in each miScript miRNA PCR Array

Control	Purpose
<i>C. elegans</i> miR-39 miScript Primer Assay	Alternative data normalization using exogenously spiked Syn-cel-miR-39 miScript miRNA Mimic
3 mature miRNA miScript Primer Assays (miR-16, miR-21, miR-191)*	Ubiquitously expressed mature miRNAs
6 snoRNA/snRNA miScript PCR Controls: SNORD61 assay SNORD95 assay SNORD96A assay SNORD68 assay† SNORD72 assay† RNU6B/RNU6-2 assay†	Data normalization using the $\Delta\Delta C_T$ method of relative quantification
miRNA reverse transcription control (miRTC miScript Primer Assay)	Assessment of reverse transcription performance
Positive PCR control (PPC)	Assessment of PCR performance

* Only included on miScript miRNA QC PCR Array.

† Only included on Pathway-Focused miScript miRNA PCR Array, miRNome miScript miRNA PCR Array, and miScript miRNA HC PCR Array.

The miScript Primer Assay for *C. elegans* miR-39 detects the miRNeasy Serum/Plasma Spike-In Control (cat. no. 219610; also called Syn-cel-miR-39 miScript miRNA Mimic [cat. no. MSY0000010]), which is a *C. elegans* miR-39 mimic. This mimic can be added to samples, particularly serum or plasma samples, to control for variations during the preparation of total RNA and subsequent steps. After purification, real-time RT-PCR detection of the miRNeasy Serum/Plasma Spike-In Control can be performed, and these results can then be used for normalization of real-time RT-PCR results for endogenous miRNAs in the sample. For more information on the use of the miRNeasy Serum/Plasma Spike-In Control, see the *miRNeasy Serum/Plasma Handbook*.

The miScript Primer Assays for miR-16, miR-21, and miR-191 are included in the miScript miRNA QC PCR Array since their targeted miRNAs are ubiquitously expressed in a wide range of cells, tissues, and body fluid samples (including whole blood, serum, and plasma). This is in contrast to the snoRNA/snRNA miScript PCR Controls whose intended targets are typically only expressed in cells, tissues, and whole blood samples. In addition, the sequences for these

mature miRNAs are conserved across a wide range of species, including all species currently supported by miScript miRNA PCR Arrays. These benefits make the miScript Primer Assays for miR-16, miR-21, and miR-191 ideal for sample quality assessment.

For accurate and reproducible results in miRNA quantification by real-time PCR, it is necessary to normalize the amount of target miRNA by using a suitable endogenous reference RNA. This approach is known as relative quantification. Normalization corrects for factors that could otherwise lead to inaccurate quantification. These factors include variation in quantity of input RNA, possible RNA degradation or presence of inhibitors in the RNA samples, and differences in sample handling. Normalization also allows results from different experiments and samples to be compared directly. miScript PCR Controls are primers designed to quantify a panel of 5 snoRNAs (SNORD61, SNORD68, SNORD72, SNORD95, and SNORD96A) and the snRNA RNU6B (RNU6-2). These controls take into consideration sequence homologies in human, mouse, rat, dog, and rhesus macaque so that the same controls can be used for all 4 species. In addition, these small RNAs have been verified to have relatively stable expression levels across tissues and cell types. As a result, miScript PCR Controls serve as normalization controls for relative quantification using the miScript PCR System. All the controls have amplification efficiencies close to 100%.

The miRTC miScript Primer Assay is an assay that assesses the performance of a reverse-transcription reaction using the miScript II RT Kit by detecting template synthesized from the kit's built-in miRNA reverse transcription control RNA (miRTC). This control monitors for any variables that may inhibit the reverse transcription reaction.

The positive PCR control (PPC) wells contain a predisposed artificial DNA sequence and the assay that detects it. This control monitors for any variables that may inhibit the PCR reaction.

Data analysis

Data analysis can then be conducted at QIAGEN's GeneGlobe Data Analysis Center using a software-based tool or with a spreadsheet-based tool that can be downloaded from a QIAGEN website.

Note: The GeneGlobe Data Analysis Center is a web resource for the analysis of real-time PCR or NGS data (www.qiagen.com/us/shop/genes-and-pathways/data-analysis-center-overview-page/?akamai-feo=off). To access the center, new users can register online. Once on the site, the data analysis software will be found under "Analysis." The miScript miRNA PCR Data Analysis spreadsheets can be found under "Product Resources/Performance Data" at www.qiagen.com/us/shop/pcr/primer-sets/miscript-mirna-pcr-arrays/#resources.

Note: If using a 384-well format, download the PCR Array 4x96 384-Well Conversion spreadsheet to dissect a 384-well dataset into the correct 4 sets of 96 genes for each of the 4 samples. For software-based data analysis, this file can be found at the GeneGlobe Data Analysis Center. For spreadsheet-based data analysis, this file can be found under "Product Resources/Performance Data" at www.qiagen.com/us/shop/pcr/primer-sets/miscript-mirna-pcr-arrays/#resources.

Template RNA requirements

Total RNA containing miRNA is the required starting material for miScript miRNA PCR Arrays. It is not necessary to enrich for small RNA. QIAGEN provides a range of solutions for purification of total RNA including miRNA (Table 2).

Table 2. Kits for purification of RNA including miRNA

Kit	Cat. no.	Starting material
miRNeasy Micro Kit	217084	Small amounts of cells and tissues
miRNeasy Mini Kit	217004	Animal/human tissues and cells
miRNeasy 96 Kit	217061	Animal/human tissues and cells
miRNeasy Serum/Plasma Kit	217184	Animal and human plasma and serum
miRNeasy FFPE Kit	217504	FFPE tissue samples
PAXgene® Tissue miRNA Kit	766134	Animal/human tissues that have been fixed and stabilized in PAXgene Tissue Containers
PAXgene Blood miRNA Kit	763134	Human blood that has been stabilized in PAXgene Blood RNA Tubes

Equipment and Reagents to Be Supplied by User

When working with chemicals, always wear a suitable lab coat, disposable gloves and protective goggles. For more information, consult the appropriate safety data sheets (SDS) available from the product supplier.

For reverse transcription

- Thin-walled, DNase-free, RNase-free PCR tubes (for 20 µl reactions)
- Ice
- Thermal cycler, heating blocks, or water baths (capable of reaching 95°C)
- Microcentrifuge

For quantitative, real-time PCR

- Real-time PCR cycler; the table on page 8 shows the appropriate real-time cycler for each array format

-
- Multichannel pipettor
 - Nuclease-free pipet tips and tubes

Protocol: Reverse Transcription for Quantitative, Real-Time PCR

Important points before starting

- The miScript II RT Kit includes 2 buffers: 5x miScript HiSpec Buffer and 5x miScript HiFlex Buffer. Only 5x miScript HiSpec Buffer should be used to prepare cDNA for mature miRNA profiling using Pathway-Focused or miRNome miScript miRNA PCR Arrays or miScript miRNA HC PCR Arrays.
- **IMPORTANT:** Only miScript HiSpec Buffer should be used to prepare cDNA for real-time PCR with miScript miRNA PCR Arrays.
- **IMPORTANT:** Do not use miScript HiFlex Buffer to prepare cDNA for real-time PCR with miScript miRNA PCR Arrays.
- **IMPORTANT:** If performing a reverse-transcription reaction for use with the miScript PreAMP PCR Kit, refer to the reverse-transcription protocol in the *miScript PreAMP Handbook*.
- Total RNA containing miRNA should be used as starting material. For RNA purification recommendations, see page 27. This protocol is for use with up to 2 µg quantifiable RNA or up to 9 µl RNA prepared from 100–200 µl serum or plasma using the miRNeasy Serum/Plasma Kit. If using higher RNA amounts, scale up the reaction linearly. Recommended starting amounts are shown in Table 3. If working with RNA for the first time, read Appendix B.
- If an miRNA-enriched fraction has been prepared separately from larger RNAs (e.g., using the miRNeasy Mini Kit), use one-tenth the recommended RNA input for total RNA. For example, if 125–250 ng total RNA is recommended for use with a Pathway-Focused miScript miRNA PCR Array, use 12.5–25 ng miRNA-enriched fraction.
- Set up all reactions on ice to minimize the risk of RNA degradation.

- Do not vortex template RNA or any of the components of the miScript II RT Kit.

PCR application	Assay	Buffer	Recommended RNA input*
Pathway profiling of mature miRNA	Pathway-Focused miScript miRNA PCR Arrays	5x miScript HiSpec Buffer	125–250 ng per RNA sample [†]
High-content pathway profiling of mature miRNA	miScript miRNA HC PCR Arrays	5x miScript HiSpec Buffer	250–500 ng per RNA sample [†]
Whole miRNome profiling of mature miRNA	miRNome miScript miRNA PCR Arrays	5x miScript HiSpec Buffer	250–500 ng per 384-well plate or per 4 x 96-well plates/ Rotor-Discs (the number of plates provided in a miRNome miScript miRNA PCR Array varies depending on the species of interest) ^{†§}

Table 3. Recommended RNA starting amounts and buffers for reverse-transcription reactions

* If the RNA sample is not limiting, use the upper amount of the recommended range.

[†] If RNA was prepared from 100–200 µl serum or plasma using the miRNeasy Serum/Plasma Kit, we recommend using 1.5 µl RNA prep (approximately one-tenth final eluate) per miScript miRNA PCR Array.

[†] If RNA was prepared from 100–200 µl serum or plasma using the miRNeasy Serum/Plasma Kit, we recommend using 1.5 µl RNA prep (approximately one-tenth final eluate) per 384-well plate or per 4 x 96-well plates/Rotor-Discs.

[§] Depending on the RNA starting amount, a single reverse-transcription reaction can provide sufficient cDNA for 8 x 384-well plates or 32 x 96-well plates/Rotor-Discs.

Procedure

1. Thaw template RNA on ice. Thaw RNase-free water, 10x miScript Nucleics Mix and 5x miScript HiSpec Buffer at room temperature (15–25°C).

Mix each solution by flicking the tubes. Centrifuge briefly to collect residual liquid from the sides of the tubes and then store on ice.

2. Prepare the reverse-transcription reaction on ice according to Table 4.

Gently mix and then store on ice. The reverse-transcription master mix contains all components required for first-strand cDNA synthesis except template RNA.

Note: miScript Reverse Transcriptase Mix should be removed from the –20°C freezer just before preparation of the master mix, gently mixed and placed on ice. It should be returned to the freezer immediately after use.

Note: If setting up more than 1 reaction, prepare a volume of master mix 10% greater than that required for the total number of reactions to be performed.

Table 4. Reverse-transcription reaction components

Component	Volume/reaction
5x miScript HiSpec Buffer	4 μ l
10x miScript Nucleic Mix	2 μ l
RNase-free water	Variable
miScript Reverse Transcriptase Mix	2 μ l
Template RNA (added in step 3)	Variable (see Table 3 for recommendations) *
Total volume	20 μl

* If RNA was prepared from 100–200 μ l serum or plasma using the miRNeasy Serum/Plasma Kit, up to 9 μ l RNA prep can be added to the reverse-transcription reaction (sufficient for 6 x 384-well plates or 24 x 96-well plates/Rotor-Discs).

3. Add template RNA to each tube containing reverse-transcription master mix. Gently mix, briefly centrifuge and then store on ice.
4. Incubate for 60 minutes at 37°C.
5. Incubate for 5 minutes at 95°C to inactivate miScript Reverse Transcriptase Mix and place on ice.
6. Dilute the cDNA in RNase-free water according to Table 5, and proceed with real-time PCR immediately.

If you wish to store the reverse-transcription reactions prior to real-time PCR, transfer the undiluted cDNA to a –20°C freezer, or dispense the diluted cDNA into 110 μ l aliquots and transfer them to a –20°C freezer.

Table 5. cDNA dilution prior to PCR

PCR application	Array	Reaction dilution
Pathway profiling	Pathway-Focused miScript miRNA PCR Arrays	Add 200 μ l RNase-free water to each 20 μ l reverse-transcription reaction
High-content pathway profiling	miScript miRNA HC PCR Arrays	Add 90 μ l RNase-free water to each 20 μ l reverse-transcription reaction
Whole miRNome profiling	miRNome miScript miRNA PCR Arrays	Dilution depends on the number of plates/Rotor-Discs in the miRNome miScript miRNA PCR Array: For 1 x 384-well plate or 4 x 96-well plates/Rotor-Discs: add 90 μ l RNase-free water to the 20 μ l reverse-transcription reaction. For 2 x 384-well plate or 8 x 96-well plates/Rotor-Discs: add 200 μ l RNase-free water to the 20 μ l reverse-transcription reaction. For 3 x 384-well plate or 12 x 96-well plates/Rotor-Discs: add 310 μ l RNase-free water to the 20 μ l reverse-transcription reaction. For 4 x 384-well plate or 16 x 96-well plates/Rotor-Discs: add 420 μ l RNase-free water to the 20 μ l reverse-transcription reaction.

Protocol: Real-Time PCR for Mature miRNA Expression Profiling

cDNA prepared using the miScript II RT Kit with miScript HiSpec Buffer (with optional preamplification, if required, using the miScript PreAMP PCR Kit) is the appropriate starting material for this protocol. This protocol enables real-time PCR profiling of mature miRNA using miScript miRNA PCR Arrays in combination with the miScript SYBR Green PCR Kit, which contains the miScript Universal Primer (reverse primer) and QuantiTect SYBR Green PCR Master Mix.

Important points before starting

- The PCR must start with an initial incubation step of 15 minutes at 95°C to activate HotStarTaq DNA Polymerase (included in 2x QuantiTect SYBR Green PCR Master Mix).
- **IMPORTANT:** Only miScript HiSpec Buffer should be used to prepare cDNA for real-time PCR with miScript miRNA PCR Arrays.
- **IMPORTANT:** Do not use miScript HiFlex Buffer to prepare cDNA for real-time PCR with miScript miRNA PCR Arrays.
- Ensure that the 20 µl cDNA synthesis reaction has been diluted appropriately. See Table 5 for recommendations. If cDNA has been preamplified using the miScript PreAMP PCR Kit with a miScript PreAMP Primer Mix, see the *miScript PreAMP Handbook* for dilution recommendations.
- Do not vortex template cDNA or any of the components of the miScript SYBR Green PCR Kit.
- The standard miScript miRNA PCR Array reaction volumes are 10 µl per well for a 384-well plate, 25 µl per well for a 96-well plate, and 20 µl per well for a 100-well Rotor-Disc.

- The miScript SYBR Green PCR Kit (200) provides sufficient reagents for 4 x 96-well arrays (110 µl residual), 2 x 384-well pathway arrays (1210 µl residual), 2 x 384-well HC/miRNome arrays (1510 µl residual), or 5 Rotor-Disc 100 arrays (110 µl residual). The miScript SYBR Green PCR Kit (1000) provides sufficient reagents for 20 x 96-well arrays (0 µl residual), 12 x 384-well pathway arrays (1100 µl residual), 13 x 384-well HC/miRNome arrays (850 µl residual), or 25 Rotor-Disc 100 arrays (0 µl residual).
- If using the iCycler iQ, iQ5, or MyiQ, well factors must be collected at the beginning of each experiment. Well factors are used to compensate for any system or pipetting nonuniformity. For details, refer to the user manual supplied with the instrument or *Technical Information: Using QuantiTect SYBR Green Kits on Bio-Rad cyclers* available at www.qiagen.com/knowledgehub.
- The miScript miRNA QC PCR Array can be used to assess the quality of cDNA samples prior to running a miScript miRNA PCR Array.

Procedure

1. Thaw 2x QuantiTect SYBR Green PCR Master Mix, 10x miScript Universal Primer, template cDNA, and RNase-free water at room temperature (15–25°C). Mix the individual solutions.
2. Prepare a reaction mix according to either Table 6 (for a Pathway-Focused miScript miRNA PCR Array) or Table 7 (for a miRNome miScript miRNA PCR Array). Mix thoroughly but gently.

Due to the hot start, it is not necessary to keep samples on ice during reaction setup or while programming the real-time cycler.

Table 6. Reaction mix for Pathway-Focused miScript miRNA PCR Arrays

Array format: Component	384-well (4 x 96) Formats E, G*	96-well Formats A, C, D, F	Rotor-Disc 100 Format R
2x QuantiTect SYBR Green PCR Master Mix [†]	550 µl	1375 µl	1100 µl
10x miScript Universal Primer	110 µl	275 µl	220 µl
RNase-free water	340 µl	1000 µl	780 µl
Template cDNA [‡]	100 µl	100 µl	100 µl
Total volume	1100 µl	2750 µl	2200 µl

* Volumes shown are sufficient for 1 cDNA template. In total, 4 cDNA templates can be analyzed on one 384-well Pathway-Focused miScript miRNA PCR Array because assays are arrayed in quadruplicate (see Figure 4).

[†] No optimization of the Mg²⁺ concentration is required. The final Mg²⁺ concentration provided by 2x QuantiTect SYBR Green PCR Master Mix gives optimal results.

[‡] Provides 0.5–1 ng cDNA per well.

Table 7. Reaction mix for miRNome miScript miRNA PCR Arrays* and miScript miRNA HC PCR Arrays[†]

Array format: Component	384-well (4 x 96) Formats E, G*	96-well Formats A, C, D, F	Rotor-Disc 100 Format R
2x QuantiTect SYBR Green PCR Master Mix [†]	2050 µl	1375 µl	1100 µl
10x miScript Universal Primer	410 µl	275 µl	220 µl
RNase-free water	1540 µl	1075 µl	855 µl
Template cDNA [§]	100 µl	25 µl	25 µl
Total volume	4100 µl	2750 µl	2200 µl

* Volumes are for a single plate or Rotor-Disc associated with a miRNome set. The number of plates in a miRNome set vary depending on the species. Scale up volumes according to the number of plates/Rotor-Discs to be run. If the miRNome set contains a plate/Rotor-Disc that is less than half full, scale down volumes accordingly.

[†] For miScript miRNA HC PCR Arrays, use the volumes shown for 384-well, Formats E, G.

[‡] No optimization of the Mg²⁺ concentration is required. The final Mg²⁺ concentration provided by 2x QuantiTect SYBR Green PCR Master Mix gives optimal results.

[§] Provides 0.5–1 ng cDNA per well.

3. Carefully remove the miScript miRNA PCR Array from its sealed bag.

Optional for 96-well and 384-well array formats: If the reaction mix is in a tube, transfer to a loading reservoir, such as the RT² PCR Array Loading Reservoir (cat. no. 338162).

4. Add reaction mix to each well of the miScript miRNA PCR Array as follows:

Note: For 384-well and 96-well array formats, a multichannel pipettor can be used to add reaction mix to the array. For the Rotor-Disc 100 format, a repeater pipettor or a QIAgility® can be used to load the array.

For 384-well miScript miRNA PCR Array: add 10 µl per well.

For 96-well miScript miRNA PCR Array: add 25 µl per well.


For Rotor-Disc miScript miRNA PCR Array: add 20 µl per well.

Loading 384-well (4 x 96) Pathway-Focused miScript miRNA PCR Arrays:


Note: Each Pathway-Focused miScript miRNA PCR Array contains 4 replicates of 96 assays that can be used for analysis of 4 samples. The spacing between the tips of standard multichannel pipettors allows rows or columns to be skipped when adding each sample. Be sure to load each sample into the correct set of wells using a multichannel pipettor and the 384EZLoad Covers (provided). Use Figure 10 as a guide. Do not reuse 384EZLoad Covers.

- 4a. Place 384EZLoad Cover 1 (white) on the plate. Add 10 µl reaction mix for sample 1 to the open wells (odd number wells of rows A, C, E, G, I, K, M, and O). Remove and discard 384EZLoad Cover 1.
- 4b. Place 384EZLoad Cover 2 (yellow) on the plate. Add 10 µl reaction mix for sample 2 to the open wells (even number wells of rows A, C, E, G, I, K, M, and O). Remove and discard 384EZLoad Cover 2.
- 4c. Place 384EZLoad Cover 3 (black) on the plate. Add 10 µl reaction mix for sample 3 to the open wells (odd number wells of rows B, D, F, H, J, L, N, and P). Remove and discard 384EZLoad Cover 3.
- 4d. Place 384EZLoad Cover 4 (red) on the plate. Add 10 µl reaction mix for sample 4 to the open wells (even number wells of rows B, D, F, H, J, L, N, and P). Remove and discard 384EZLoad Cover 4.


Cover 1 (white) for sample 1

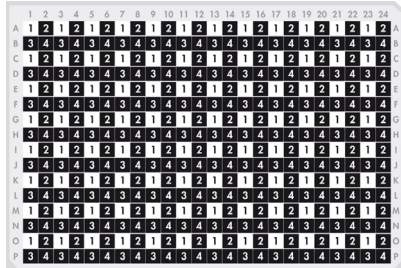

Cover 2 (yellow) for sample 2

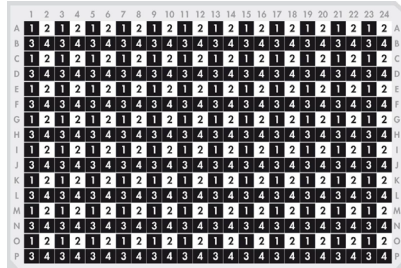

Cover 3 (black) for sample 3

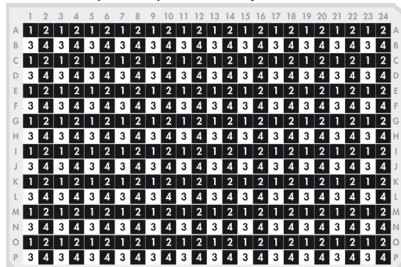

Cover 4 (red) for sample 4

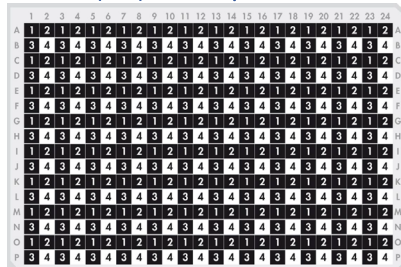


Figure 10. Loading Pathway-Focused miScript miRNA PCR Arrays, plate format E, G (384-well).

Add 10 µl reaction mix for each of 4 samples into the staggered wells with the same number as indicated in the figure.

- Carefully, tightly seal the miScript miRNA PCR Array with Optical Thin-Wall 8-Cap Strips (Formats A and D), Optical Adhesive Film (Formats C, E, F, and G), or Rotor-Disc Heat-Sealing Film (Format R).
- Centrifuge the PCR plate for 1 minute at 1000 g at room temperature (15–25°C) to remove bubbles.

Note: This step is not necessary for reactions set up in Rotor-Discs.
- Program the real-time cyclers according to Table 8.

Note: Perform dissociation curve analysis of the PCR product(s) to verify their specificity and identity. Dissociation curve analysis is an analysis step built into the software of real-time cyclers. Follow the instructions provided by the supplier.

Table 8. Cycling conditions for real-time PCR

Step	Time	Temperature	Additional comments
PCR Initial activation step	15 min	95°C	HotStarTaq DNA Polymerase is activated by this heating step.
3-step cycling:^{**†}			
Denaturation	15 s	94°C	
Annealing	30 s	55°C	
Extension [§]	30 s	70°C	Perform fluorescence data collection.
Cycle number	40 cycles		Cycle number depends on the amount of template cDNA and abundance of the target.

* For Bio-Rad models CFX96 and CFX384: adjust the ramp rate to 1°C/s.

† For Eppendorf Mastercycler ep realplex models 2, 2S, 4, and 4S: for the Silver Thermoblock, adjust the ramp rate to 26%; for the Aluminum Thermoblock, adjust the ramp rate to 35%.

‡ If using a Roche LightCycler 480, adjust the ramp rate to 1°C/s.

§ Due to software requirements, the fluorescence detection step must be at least 30 s with the ABI PRISM[®] 7000 or 34 s with the Applied Biosystems 7300 and 7500.

¶ If using a Roche LightCycler 480, use 45 cycles.

8. Place the plate/Rotor-Disc in the real-time cycler and start the cycling program.

9. Perform data analysis.

Protocol: Data Analysis

This protocol describes the steps for analysis of data from miScript miRNA PCR Arrays. The first steps should be performed by the user. The later steps can be performed at QIAGEN's GeneGlobe Data Analysis Center using a software-based tool or with a spreadsheet-based tool that can be downloaded from a QIAGEN website.

Note: The GeneGlobe Data Analysis Center is a web resource for the analysis of real-time PCR or NGS data (www.qiagen.com/us/shop/genes-and-pathways/data-analysis-center-overview-page/?akamai-feo=off). To access the center, new users can register online. Once on the site, the data analysis software will be found under "Analysis." The miScript miRNA PCR Data Analysis spreadsheet can be found under "Product Resources/Performance Data" at www.qiagen.com/us/shop/pcr/primer-sets/miscript-mirna-pcr-arrays/#resources.

Note: If using a 384-well format, download the PCR Array 4x96 384-Well Conversion spreadsheet to dissect a 384-well dataset into the correct 4 sets of 96 genes for each of the 4 samples. For software-based data analysis, this file can be found at the GeneGlobe Data Analysis Center. For spreadsheet-based data analysis, this file can be found under "Product Resources/Performance Data" at www.qiagen.com/us/shop/pcr/primer-sets/miscript-mirna-pcr-arrays/#resources.

Important points before starting

- Text marked with a ■ denotes instructions for 96-well and 384-well plates (formats A, C, D, E, F, and G); text marked with a ▲ denotes instructions for 100-well Rotor-Discs (format R).

Procedure

Steps performed by the user

1. Define the baseline.

The baseline is the noise level in early cycles, where there is no detectable increase in fluorescence due to PCR products.

■ Use the “Linear View” of the amplification plot to determine the earliest visible amplification. Set the baseline from cycle 2 to 2 cycles before the earliest visible amplification. Do not use greater than cycle 15. The number of cycles used to calculate the baseline can be changed and should be reduced if high template amounts are used. For more information regarding real-time PCR data output, refer to Appendix A.

▲ For the Rotor-Gene Q, we recommend using the “Dynamic Tube” setting along with the “Slope Correct” and/or “Ignore First” settings. For more information, refer to the Rotor-Gene Q User Manual.

Note: Ensure that baseline settings are the same across all PCR runs associated with the same experiment to allow comparison of results.

2. Define the threshold.

The threshold should be set using a logarithmic amplification plot so that the log-linear range of the curve can be easily identified. Using the “Log View” of the amplification plot, place the threshold above the background signal but within the lower half of the log-linear range of the amplification plot. The threshold should never be set in the plateau phase. The absolute position of the threshold is less critical than its consistent position across PCR runs.

■ Various PCR instruments (such as Applied Biosystems models 7500 and ViiA 7, and Stratagene models Mx3005P and Mx3000P) may require adjustment of the default “Manual C_T” threshold value of 0.2 to a lower value in order to analyze the data properly. Use a value of 0.02 as a starting point.

▲ For the Rotor-Gene Q, we recommend a C_T threshold value of approximately 0.02 in order to analyze the data properly.

Note: Ensure that threshold settings are the same across all PCR runs in the same analysis to allow comparison of results.

3. Export C_T values according to the manual supplied with the real-time PCR instrument.

Protocol: cDNA Quality Control Prior to Profiling Mature miRNA

Use of the miScript miRNA QC PCR Array enables testing of the quality of cDNA prepared using miScript HiSpec Buffer, saving time and reagents. Only 5 µl diluted cDNA sample is required for quality control using the miScript miRNA QC PCR Array. cDNA prepared using the miScript II RT Kit with miScript HiSpec Buffer is the appropriate starting material for this protocol. This protocol describes quality control of multiple cDNA samples using the miScript SYBR Green PCR Kit and miScript miRNA QC PCR Array prior to miRNA profiling using miScript miRNA PCR Arrays. In total, 32 cDNA samples can be analyzed on one 384-well miScript miRNA QC PCR Array, 8 cDNA samples can be analyzed on one 96-well miScript miRNA QC PCR Array, and 8 cDNA samples can be analyzed on one Rotor-Disc 100 miScript miRNA QC PCR Array.

Important points before starting

- The PCR must start with an initial incubation step of 15 minutes at 95°C to activate HotStarTaq DNA Polymerase (included in 2x QuantiTect SYBR Green PCR Master Mix).
- Only cDNA template prepared using miScript HiSpec Buffer should be used with this protocol.
- Ensure cDNA samples were prepared according to the protocol on page 29, and that the 20 µl cDNA synthesis reaction has been diluted appropriately (see Table 5).
- Do not vortex template cDNA or any of the components of the miScript SYBR Green PCR Kit.
- The standard miScript miRNA PCR Array reaction volumes are 10 µl per well for a 384-well plate, 25 µl per well for a 96-well plate, and 20 µl per well for a 100-well Rotor-Disc.

- If using the iCycler iQ, iQ5, or MyiQ, well factors must be collected at the beginning of each experiment. Well factors are used to compensate for any system or pipetting nonuniformity. For details, refer to the user manual supplied with the instrument or *Technical Information: Using QuantiTect SYBR Green Kits on Bio-Rad cyclers* available at www.qiagen.com/knowledgehub.

Procedure

1. Thaw 2x QuantiTect SYBR Green PCR Master Mix, 10x miScript Universal Primer, template cDNA, and RNase-free water at room temperature (15–25°C). Mix the individual solutions.
2. Prepare a reaction mix according to Table 9. Mix gently and thoroughly.

Due to the hot start, it is not necessary to keep samples on ice during the reaction setup or while programming the real-time cycler.

Table 9. Reaction mix for miScript miRNA QC PCR Arrays*

Array format: Component	384-well (4 x 96) Formats E, G†	96-well Formats A, C, D, F†	Rotor-Disc 100 Format R†
2x QuantiTect SYBR Green PCR Master Mix‡	75 µl	175 µl	150 µl
10x miScript Universal Primer	15 µl	35 µl	30 µl
RNase-free water	55 µl	135 µl	115 µl
Template cDNA	5 µl	5 µl	5 µl
Total volume	150 µl	350 µl	300 µl

* These volumes provide a 10 µl per well reaction volume for a 384-well plate, 25 µl per well reaction volume for a 96-well plate, and 20 µl per well reaction volume for a 100-well Rotor-Disc.

† Volumes shown are sufficient for 1 cDNA template. In total, 8 cDNA samples can be analyzed on one 96-well miScript miRNA QC PCR Array, 8 cDNA samples can be analyzed on one Rotor-Disc 100 miScript miRNA QC PCR Array, and 32 cDNA samples can be analyzed on one 384-well miScript miRNA QC PCR Array.

‡ No optimization of the Mg²⁺ concentration is required. The final Mg²⁺ concentration provided by 2x QuantiTect SYBR Green PCR Master Mix gives optimal results.

3. Carefully remove the miScript miRNA QC PCR Array from its sealed bag.
4. Add reaction mix to the wells of the miScript miRNA QC PCR Array as follows:

For 384-well miScript miRNA PCR Array: add 10 µl per well.

For 96-well miScript miRNA PCR Array: add 25 µl per well.

For Rotor-Disc 100 miScript miRNA PCR Array: add 20 µl per well.

- Carefully, tightly seal the miScript miRNA QC PCR Array with Optical Thin-Wall 8-Cap Strips (Formats A and D), Optical Adhesive Film (Formats C, E, F, and G), or Rotor-Disc Heat-Sealing Film (Format R).
- Centrifuge the PCR plate for 1 minute at 1000 g at room temperature (15–25°C) to remove bubbles.

Note: This step is not necessary for reactions set up in Rotor-Discs.

- Program the real-time cycler according to Table 10.

Note: Perform dissociation curve analysis of the PCR product(s) to verify their specificity and identity. Dissociation curve analysis is an analysis step built into the software of real-time cyclers. Follow the instructions provided by the supplier.

Table 10. Cycling conditions for real-time PCR

Step	Time	Temperature	Additional comments
PCR Initial activation step	15 min	95°C	HotStarTaq DNA Polymerase is activated by this heating step.
3-step cycling:**			
Denaturation	15 s	94°C	
Annealing	30 s	55°C	
Extension [§]	30 s	70°C	Perform fluorescence data collection.
Cycle number	40 cycles		Cycle number depends on the amount of template cDNA and abundance of the target.

* For Bio-Rad models CFX96 and CFX384: adjust the ramp rate to 1°C/s.

† For Eppendorf Mastercycler ep realplex models 2, 2S, 4, and 4S: for the Silver Thermoblock, adjust the ramp rate to 26%; for the Aluminum Thermoblock, adjust the ramp rate to 35%.

‡ If using a Roche LightCycler 480, adjust the ramp rate to 1°C/s.

§ Due to software requirements, the fluorescence detection step must be at least 30 s with the ABI PRISM 7000 or 34 s with the Applied Biosystems 7300 and 7500.

¶ If using a Roche LightCycler 480, use 45 cycles.

7. Place the plate/Rotor-Disc in the real-time cycler and start the cycling program.
8. Perform data analysis.

Protocol: Data Analysis for Quality Control Using miScript miRNA QC PCR Arrays

This protocol describes the steps for analysis of data from miScript miRNA QC PCR Arrays. The first steps should be performed by the user. The later steps are performed at QIAGEN's GeneGlobe Data Analysis Center using a software-based tool or with a spreadsheet-based tool that can be downloaded from a QIAGEN website.

Important point before starting

- Text marked with a ■ denotes instructions for 96-well and 384-well plates (formats A, C, D, E, F, and G); text marked with a ▲ denotes instructions for 100-well Rotor-Discs (format R).

Procedure

Steps performed by the user

1. Define the baseline.

■ Use the "Linear View" of the amplification plot to determine the earliest visible amplification. Set the baseline from cycle 2 to 2 cycles before the earliest visible amplification. Do not use greater than cycle 15. The number of cycles used to calculate the baseline can be changed and should be reduced if high template amounts are used. For more information regarding real-time PCR data output, refer to Appendix A.

▲ For the Rotor-Gene Q, we recommend using the "Dynamic Tube" setting along with the "Slope Correct" and/or "Ignore First" settings. For more information, refer to the *Rotor-Gene Q User Manual*.

Note: Ensure that baseline settings are the same across all PCR runs in the same analysis to allow comparison of results.

2. Define the threshold.

The threshold should be set using a logarithmic amplification plot so that the log-linear range of the curve can be easily identified. Using the “Log View” of the amplification plot, place the threshold above the background signal but within the lower half of the log-linear range of the amplification plot. The threshold should never be set in the plateau phase. The absolute position of the threshold is less critical than its consistent position across PCR runs.

■ Various PCR instruments (such as Applied Biosystems models 7500 and ViiA 7, and Stratagene models Mx3005P and Mx3000P) may require adjustment of the default “Manual C_T ” threshold value of 0.2 to a lower value in order to analyze the data properly. Use a value of 0.02 as a starting point.

▲ For the Rotor-Gene Q, we recommend a C_T threshold value of approximately 0.02 in order to analyze the data properly.

Note: Ensure that threshold settings are the same across all PCR runs in the same analysis to allow comparison of results.

3. Export C_T values according to the manual supplied with the real-time PCR instrument.
4. Data analysis can then be conducted at QIAGEN’s GeneGlobe Data Analysis Center using a software-based tool or with a spreadsheet-based tool that can be downloaded from a QIAGEN website.

Note: The GeneGlobe Data Analysis Center is a web resource for the analysis of real-time PCR or NGS data (www.qiagen.com/us/shop/genes-and-pathways/data-analysis-center-overview-page/?akamai-feo=off). To access the center, new users can register online. Once on the site, the data analysis software will be found under “Analysis.” The miScript miRNA QC PCR Data Analysis spreadsheet can be found under

“Product Resources/Performance Data” at www.qiagen.com/us/shop/pcr/primer-sets/miscript-mirna-pcr-arrays/#resources.

Note: If using a 384-well format, download the PCR Array 4x96 384-Well Conversion spreadsheet to dissect a 384-well dataset into the correct 4 sets of 96 genes for each of the 4 samples. For software-based data analysis, this file can be found at the GeneGlobe Data Analysis Center. For spreadsheet-based data analysis, this file can be found under “Product Resources/Performance Data” at www.qiagen.com/us/shop/pcr/primer-sets/miscript-mirna-pcr-arrays/#resources.

Troubleshooting Guide

This troubleshooting guide may be helpful in solving any problems that may arise. For more information, see also the Frequently Asked Questions page at our Technical Support Center: www.qiagen.com/us/shop/pcr/primer-sets/miscript-mirna-pcr-arrays/#resources. The scientists in QIAGEN Technical Services are always happy to answer any questions you may have about either the information and protocols in this handbook or sample and assay technologies (for contact information, see back cover or visit www.qiagen.com).

Comments and suggestions

Evidence of poor reverse transcription efficiency (value of $AVG C_T^{miRQC} - AVG C_T^{PPC} > 7$)

- | | | |
|----|---|--|
| a) | Poor quality RNA | Check the $A_{260}:A_{280}$ and $A_{260}:A_{230}$ ratios of the RNA samples. Be sure to perform the dilutions for spectrophotometry in RNase-free 10 mM Tris-Cl, pH 7.5. If necessary, repurify RNA with a spin-column based clean up method, such as the miRNeasy Mini Kit (cat. no. 217004). |
| b) | Calculation did not include correction factor | If using miRNome miScript miRNA PCR Arrays, miScript miRNA HC PCR Arrays, or miScript miRNA QC PCR Arrays, be sure to include the correction factors detailed in Table 9 or Table 10. |

Evidence of poor, overall PCR amplification efficiency ($AVG C_T^{PPC}$ varies by more than 2 across arrays and/or is greater than 21 for 96-well and 384-well plates or 17 for 100-well Rotor-Discs)

- | | | |
|----|--|---|
| a) | Variation in instrument sensitivity | Different instruments have different levels of sensitivity. If an average C_T^{PPC} value of 19 ± 2 for 96-well and 384-well plates or 15 ± 2 for 100-well Rotor-Discs is difficult to obtain for the instrument used, the observed average C_T^{PPC} value should be acceptable as long as it does not vary by more than 2 cycles between arrays being compared. |
| b) | HotStarTaq DNA Polymerase not activated with a hot start | Be sure that the initial heat activation step at 95°C took place for 15 minutes, and that all other cycle parameters were performed according to the protocol. |
| c) | Poor quality RNA that may contain PCR inhibitors | Check the $A_{260}:A_{280}$ and $A_{260}:A_{230}$ ratios of the RNA samples. Be sure to perform the dilutions for spectrophotometry in RNase-free 10 mM Tris-Cl, pH 7.5. If necessary, repurify RNA with a spin-column based clean up method, such as the miRNeasy Mini Kit (cat. no. 217004). |

Comments and suggestions

No product, or product detected late in real-time PCR (indicative of problems occurring during reverse transcription)

- | | | |
|----|---|---|
| a) | Pipetting error or missing reagent when setting up reverse-transcription reaction | Check the pipets used for experimental setup. Mix all reagents well after thawing and repeat the reverse-transcription reaction. |
| b) | Incorrect setup of reverse-transcription reaction | Be sure to set up the reaction on ice. |
| c) | Poor quality or incorrect amount of template RNA for reverse-transcription reaction | Check the concentration, integrity, and purity of the template RNA before starting the protocol. Mix well after thawing the template RNA. Even minute amounts of RNases can affect synthesis of cDNA and sensitivity in RT-PCR, particularly with small amounts of RNA. |
| d) | RNA concentration too high or too low | See Table 3 for recommended RNA amounts. |
| e) | RNA denatured | Denaturation of the template RNA is not necessary. If denaturation was performed, the integrity of the RNA may be affected. |
| f) | Incubation temperature too high | Reverse transcription should be carried out at 37°C. Higher temperatures may reduce the length of cDNA products or the activity of miScript Reverse Transcriptase Mix. Check the temperature of your heating block or water bath. |

No product, or product detected late in real-time PCR, or only primer-dimers detected (indicative of problems occurring during real-time PCR)

- | | | |
|----|--|--|
| a) | PCR annealing time too short | Use the annealing time specified in the protocol. |
| b) | PCR extension time too short | Use the extension time specified in the protocol. |
| c) | Pipetting error or missing reagent when setting up PCR | Check the concentrations and storage conditions of reagents, including primers and cDNA. |
| d) | HotStarTaq DNA Polymerase not activated with a hot start | Ensure that the cycling program includes the hot start activation step for HotStarTaq DNA polymerase; for details, check the protocol. |
| e) | No detection activated | Check that fluorescence detection was activated in the cycling program. |

Comments and suggestions

- | | | |
|----|--------------------------------|--|
| f) | Wrong detection step | Ensure that fluorescence detection takes place during the extension step of the PCR cycling program. |
| g) | Wrong dye layer/filter chosen | Ensure that the appropriate layer/filter is activated. |
| h) | Insufficient starting template | Increase the amount of template cDNA. |

No linearity in ratio of C_r value/crossing point to log of the template amount

- | | | |
|----|--------------------------|--|
| a) | Template amount too high | Do not exceed maximum recommended amounts of template cDNA. For details, see the protocol. |
| b) | Template amount too low | Increase amount of template RNA. |

Varying fluorescence intensity

- | | | |
|----|--|---|
| a) | Real-time cyclers contaminated | Decontaminate the real-time cyclers according to the supplier's instructions. |
| b) | Real-time cyclers no longer calibrated | Recalibrate the real-time cyclers according to the supplier's instructions. |

Appendix A: Real-Time PCR Data Output and Dissociation Curve Analysis

In a typical amplification plot resulting from a real-time PCR reaction, fluorescence is plotted against the number of cycles, producing sigmoidal-shaped plots (when using a linear scale). The C_T serves as a tool for calculation of the starting template amount in each sample. This is the cycle in which there is the first detectable increase in fluorescence. There may be variation in how determination of C_T values is carried out depending on the real-time PCR cycler that is used.

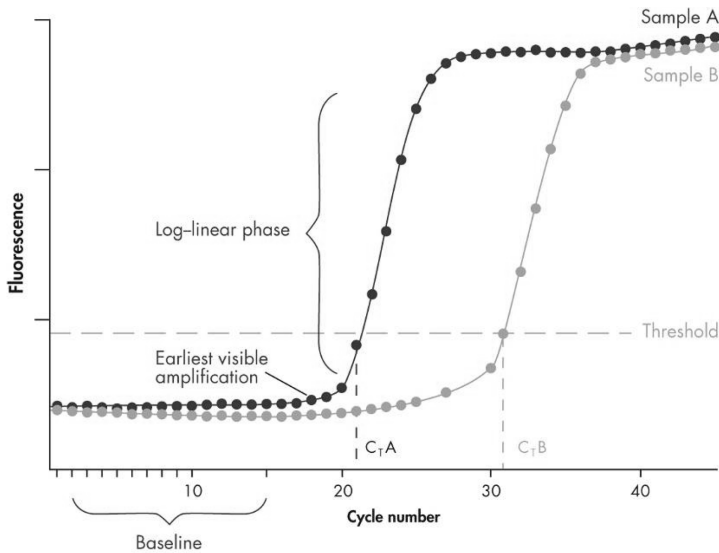


Figure 11. Amplification plot. Amplification plots showing increases in fluorescence from 2 samples (A and B). Sample A contains a higher amount of starting template than sample B.

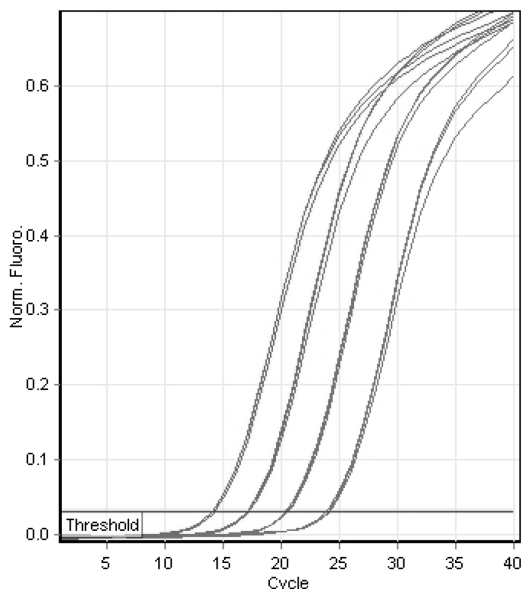


Figure 12. Typical amplification plot. Amplification plot after quantification of a range of amounts of miR-21. Real-time PCR was performed using the Rotor-Gene Q.

Dissociation curve analysis

A dissociation curve analysis of PCR product(s) may be optionally performed to aid in verifying their specificity and identity. Dissociation curve analysis is an analysis step built into the software of real-time cyclers. Follow instructions provided by the supplier.

To carry out dissociation curve analysis, the temperature is increased very slowly from a low temperature (e.g., 65°C) to a high temperature (e.g., 95°C). At low temperatures, PCR products are double stranded, so SYBR Green I dye binds to them and fluorescence is high. However at high temperatures, PCR products are denatured, resulting in rapid decreases in fluorescence.

The fluorescence is measured continuously as the temperature is increased and the fluorescence values are plotted against temperature. A curve is produced, because fluorescence decreases slightly through the lower end of the temperature range, but decreases much more rapidly at higher temperatures as the dissociation temperatures of nonspecific and specific PCR products are reached. The detection systems calculate the first derivatives of the curves, resulting in curves with peaks at the respective T_m s (Figures 13 and 14). Curves with peaks at a T_m lower than that of the specific PCR product indicate the formation of primer-dimers, while diverse peaks with different T_m s or plateaus indicate production of nonspecific products or a smear.

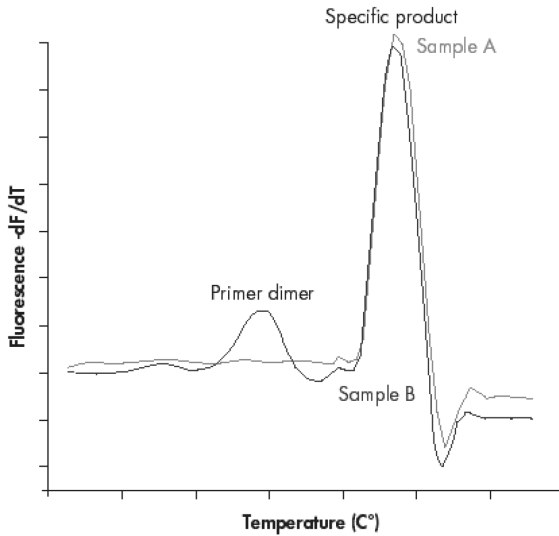


Figure 13. Dissociation curve analysis. Dissociation curve analysis of 2 samples (A and B). Sample A yields only 1 peak, resulting from the specific amplification product (primer-dimers not coamplified). Sample B shows a peak from the specific product and a peak at a lower temperature from amplification of primer-dimers.

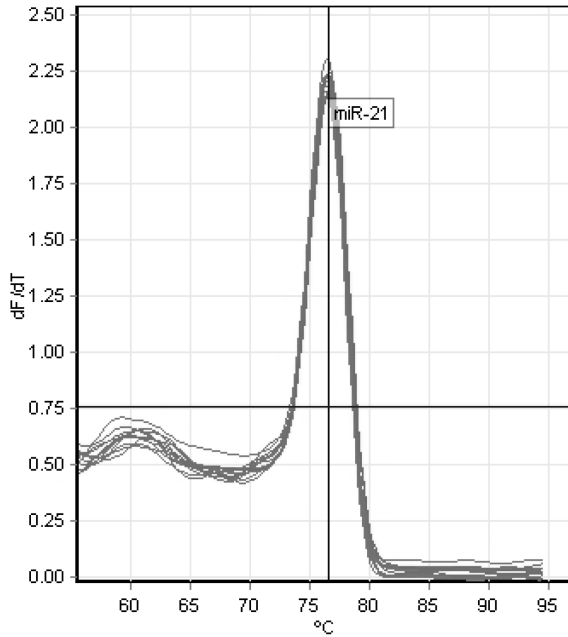


Figure 14. miRNA dissociation curve. Dissociation curve analysis of an miR-21 PCR product showing a single peak from the specific amplification product. Dissociation curve analysis was performed using the Rotor-Gene Q.

Appendix B: General Remarks on Handling RNA

Handling RNA

Ribonucleases (RNases) are very stable and active enzymes that generally do not require cofactors to function. Since RNases are difficult to inactivate and even minute amounts are sufficient to degrade RNA, do not use any plasticware or glassware without first eliminating possible RNase contamination. Care should be taken to avoid inadvertently introducing RNases into the RNA sample during or after the purification procedure. In order to create and maintain an RNase-free environment, the following precautions must be taken during pretreatment and use of disposable and nondisposable vessels and solutions while working with RNA.

General handling

Proper microbiological, aseptic technique should always be used when working with RNA. Hands and dust particles may carry bacteria and molds and are the most common sources of RNase contamination. Always wear latex or vinyl gloves while handling reagents and RNA samples to prevent RNase contamination from the surface of the skin or from dusty laboratory equipment. Change gloves frequently and keep tubes closed whenever possible. Keep purified RNA on ice when aliquots are pipetted for downstream applications.

To remove RNase contamination from bench surfaces, use of nondisposable plasticware, laboratory equipment (e.g., pipets and electrophoresis tanks) is recommended. RNase contamination can alternatively be removed using general laboratory reagents. To decontaminate plasticware, rinse with 0.1 M NaOH, 1 mM EDTA* followed by RNase-free water (see "Solutions"), or rinse with chloroform* if the plasticware is chloroform-resistant. To

* When working with chemicals, always wear a suitable lab coat, disposable gloves and protective goggles. For more information, consult the appropriate material data sheets (MSDSs), available from the product supplier.

decontaminate electrophoresis tanks, clean with detergent (e.g., 0.5% SDS),* rinse with RNase-free water, rinse with ethanol (if the tanks are ethanol-resistant), and allow to dry.

Disposable plasticware

The use of sterile, disposable polypropylene tubes is recommended throughout the procedure. These tubes are generally RNase-free and do not require pretreatment to inactivate RNases.

Glassware

Glassware should be treated before use to ensure that it is RNase-free. Glassware used for RNA work should be cleaned with a detergent,* thoroughly rinsed, and oven baked at 240°C for 4 hours or more (overnight, if more convenient) before use. Autoclaving alone will not fully inactivate many RNases. Alternatively, glassware can be treated with DEPC* (diethyl pyrocarbonate), as described in “Solutions” below.

Solutions

Note: QIAGEN solutions, such as miScript Nucleics Mix, miScript HiFlex Buffer, miScript HiSpec Buffer, and RNase-free water, are guaranteed RNase-free without using DEPC treatment and are therefore free of any DEPC contamination.

Solutions (water and other solutions) should be treated with 0.1% DEPC. DEPC is a strong, but not absolute, inhibitor of RNases. It is commonly used at a concentration of 0.1% to inactivate RNases on glass or plasticware or to create RNase-free solutions and water. DEPC inactivates RNases by covalent modification. Add 0.1 ml DEPC to 100 ml of the solution to be treated and shake vigorously to bring the DEPC into solution. Let the solution incubate for 12 hours at 37°C. Autoclave for 15 minutes to remove any trace of DEPC. DEPC will react with primary amines and cannot be used directly to treat Tris* buffers. DEPC is highly unstable in the

presence of Tris buffers and decomposes rapidly into ethanol and CO₂. When preparing Tris buffers, treat water with DEPC first, and then dissolve Tris to make the appropriate buffer. Trace amounts of DEPC will modify purine residues in RNA by carbethoxylation. Carbethoxylated RNA is translated with very low efficiency in cell-free systems. However, its ability to form DNA:RNA or RNA:RNA hybrids is not seriously affected unless a large fraction of the purine residues have been modified. Residual DEPC must always be eliminated from solutions or vessels by autoclaving or heating to 100°C for 15 minutes.

Appendix C: Preparation, Quantification, and Storage of RNA

RNA preparation and quality

Since PCR consists of multiple rounds of enzymatic reactions, it is more sensitive to impurities such as proteins, phenol/chloroform, salts, and EDTA than single-step enzyme-catalyzed reactions. Purity of nucleic acid templates is particularly important for real-time PCR, since contaminants can interfere with fluorescence detection. See Table 2 for kits recommended for the purification of total RNA that includes miRNA. For more information about kits for miRNA purification, visit www.qiagen.com/miRNA.

Determining concentration and purity of RNA

The concentration of RNA should be determined by measuring the absorbance at 260 nm (A_{260}) in a spectrophotometer. The sample should be diluted in water since the relationship between absorbance and concentration (A_{260} reading of 1 = 40 $\mu\text{g}/\text{ml}$ RNA) is based on an extinction coefficient calculated for RNA in water. To ensure significance, readings should fall between 0.15 and 1.0.

Note that absorbance measurements cannot discriminate between DNA and RNA. Depending on the method used for RNA preparation, RNA may be contaminated with DNA, and this will result in misleadingly high A_{260} values.

The ratios between the absorbance values at 260 nm and 280 nm and at 260 nm and 230 nm give an estimate of RNA purity. To determine RNA purity, we recommend measuring

absorbance in 10 mM Tris-Cl, pH 7.5. * Pure RNA has an A_{260}/A_{280} ratio of 1.9–2.1 and an A_{260}/A_{230} ratio of 2.0–2.2.† Lower ratios indicate the presence of contaminants such as proteins.

Storage of RNA

Purified RNA should be stored at -20°C or -70°C in RNase-free water. When purified using QIAGEN systems, no degradation is detectable for at least 1 year under these conditions. Diluted solutions of RNA (e.g., dilution series used as standards) should be stored in aliquots and thawed once only. We recommend storage of aliquots in siliconized tubes where possible. This avoids adsorption of the RNA to the tube walls, which would reduce the concentration of RNA in solution.

DNA contamination in RNA samples purified from different tissues

Depending on the type of tissue used as starting material for RNA purification, a fluorescent signal may also be generated in “no RT” control reactions. When RNA is purified from tissues that contain large amounts of DNA, such as spleen or thymus, the level of DNase treatment required may be higher than for other tissues. For such tissues, we recommend performing a DNase digestion (using the QIAGEN RNase-Free DNase Set, cat. no. 79254) when using the miRNeasy Mini and miRNeasy 96 Kits for RNA purification.

* When working with chemicals, always wear a suitable lab coat, disposable gloves, and protective goggles. For more information, consult the appropriate material safety data sheets (MSDSs), available from the product supplier.

† $A_{260}:A_{280}$ values up to 2.3 are routinely obtained for pure RNA (in 10 mM Tris-Cl, pH 7.5) with some spectrophotometers.

Ordering Information

Product	Contents	Cat. no.
miScript II RT Kit (12)	For 12 cDNA synthesis reactions: miScript Reverse Transcriptase Mix, 10x miScript Nucleics Mix, 5x miScript HiSpec Buffer, 5x miScript HiFlex Buffer, RNase-Free Water	218160
miScript II RT Kit (50)	For 50 cDNA synthesis reactions: miScript Reverse Transcriptase Mix, 10x miScript Nucleics Mix, 5x miScript HiSpec Buffer, 5x miScript HiFlex Buffer, RNase-Free Water	218161
miScript SYBR Green PCR Kit (200)	For 200 reactions: QuantiTect SYBR Green PCR Master Mix, miScript Universal Primer	218073
miScript SYBR Green PCR Kit (1000)	For 1000 reactions: QuantiTect SYBR Green PCR Master Mix, miScript Universal Primer	218075
miScript Primer Assay (100)	10x miScript Primer Assay (contains one miRNA-specific primer)	Varies*
Pathway-Focused miScript miRNA PCR Array	Array of assays for a pathway, disease, or gene family for human, mouse, rat, dog, or rhesus macaque miRNAs; available in 96-well, 384- well, or Rotor-Disc 100 format	Varies

*Visit www.qiagen.com/GeneGlobe to search for and order these products.

miScript miRNA HC PCR Array	High-content array of assays for a pathway, disease, or gene family miRNAs; available in 384-well format	Varies
miRNome miScript miRNA PCR Array	Array of assays for the complete human, mouse, rat, dog, or rhesus macaque miRNome; available in 96-well, 384-well, or Rotor-Disc 100 format	Varies
miScript miRNA QC PCR Array	Array of quality control assays for human, mouse, rat, dog, or rhesus macaque miRNAs; available in 96-well, 384-well, or Rotor-Disc 100 format	Varies
RT ² PCR Array Loading Reservoir	12 x 5 ml capacity, irradiation-sterilized loading reservoirs	338162
384EZLoad Covers	Pack of 4 color-coded covers for loading 384-well plates	338125
Related products		
miScript PreAMP PCR Kit (12)	For 12 preamplification reactions: 5x miScript PreAMP Buffer, HotStarTaq DNA Polymerase (2 U/μl), miScript PreAMP Universal Primer, 4 miScript Primer Assays, RNase-Free Water	331451
miScript PreAMP PCR Kit (60)	For 60 preamplification reactions: 5x miScript PreAMP Buffer, HotStarTaq DNA Polymerase (2 U/μl), miScript PreAMP Universal Primer, 4 miScript Primer Assays, RNase-Free Water	331452
miScript PreAMP Primer Mix	Primer mix for preamplification; for use with corresponding miScript miRNA PCR Array	Varies
miRNeasy Micro Kit (50)	For 50 total RNA preps: 50 RNeasy [®] MinElute [®] Spin Columns, Collection Tubes (1.5 ml and 2 ml), QIAzol [®] Lysis	217084

	Reagent, RNase-Free Reagents and Buffers	
miRNeasy Mini Kit (50)	For 50 preps: 50 RNeasy Mini Spin Columns, Collection Tubes (1.5 ml and 2 ml), QIAzol Lysis Reagent, RNase-Free Reagents and Buffers	217004
miRNeasy 96 Kit (4)	For 4 x 96 preps: 4 RNeasy 96 plates, Collection Microtubes (racked), Elution Microtubes CL, Caps, S-Blocks, AirPore Tape Sheets, QIAzol Lysis Reagent, RNase-Free Reagents and Buffers	217061
miRNeasy Serum/Plasma Kit (50)	For 50 total RNA preps: 50 RNeasy MinElute Spin Columns, Collection Tubes (1.5 ml and 2 ml), QIAzol Lysis Reagent, Ce_miR-39_1 miScript Primer Assay, RNase-free Reagents and Buffers	217184
miRNeasy Serum/Plasma Spike-In Control	10 pmol lyophilized <i>C. elegans</i> miR-39 miRNA mimic	219610
miRNeasy FFPE Kit (50)	50 RNeasy MinElute Spin Columns, Collection Tubes, Proteinase K, RNase-Free DNase I, DNase Booster Buffer, RNase-Free Buffers, RNase-Free Water	217504
PAXgene Tissue miRNA Kit (50)	For 50 RNA preps: PAXgene RNA MinElute Spin Columns, PAXgene Shredder Spin Columns, Processing Tubes, Microcentrifuge Tubes, Carrier RNA, RNase-Free DNase, and RNase-Free Buffers; to be used with PAXgene Tissue Containers	766134
PAXgene Tissue Containers (10)	For collection, fixation, and stabilization of 10 samples: 10 Prefilled Reagent Containers,	765112

PAXgene Blood miRNA Kit (50)

containing PAXgene Tissue Fix and PAXgene Tissue Stabilizer

For 50 RNA preps: PAXgene Spin Columns, PAXgene Shredder Spin Columns, Processing Tubes, Microcentrifuge Tubes, RNase-Free DNase, RNase-Free Reagents and Buffers; to be used with PAXgene Blood RNA Tubes (available from BD, cat. no. 762165)

763134

For up-to-date licensing information and product-specific disclaimers, see the respective QIAGEN kit handbook or user manual. QIAGEN kit handbooks and user manuals are available at www.qiagen.com or can be requested from QIAGEN Technical Services or your local distributor.

Document Revision History

R3	Updated template.
11/2018	Updated data analysis procedure.

Limited License Agreement for the miScript PCR System

Use of this product signifies the agreement of any purchaser or user of the product to the following terms:

1. The product may be used solely in accordance with the protocols provided with the product and this handbook and for use with components contained in the kit only. QIAGEN grants no license under any of its intellectual property to use or incorporate the enclosed components of this kit with any components not included within this kit except as described in the protocols provided with the product, this handbook, and additional protocols available at www.qiagen.com. Some of these additional protocols have been provided by QIAGEN users for QIAGEN users. These protocols have not been thoroughly tested or optimized by QIAGEN. QIAGEN neither guarantees them nor warrants that they do not infringe the rights of third-parties.
2. Other than expressly stated licenses, QIAGEN makes no warranty that this kit and/or its use(s) do not infringe the rights of third-parties.
3. This kit and its components are licensed for one-time use and may not be reused, refurbished, or resold.
4. QIAGEN specifically disclaims any other licenses, expressed or implied other than those expressly stated.
5. The purchaser and user of the kit agree not to take or permit anyone else to take any steps that could lead to or facilitate any acts prohibited above. QIAGEN may enforce the prohibitions of this Limited License Agreement in any Court, and shall recover all its investigative and Court costs, including attorney fees, in any action to enforce this Limited License Agreement or any of its intellectual property rights relating to the kit and/or its components.

For updated license terms, see www.qiagen.com.

Trademarks: QIAGEN®, QIAzol®, QIAgility®, HotStarTaq®, miScript®, QuantiTect®, RNeasy®, Rotor-Gene®, Rotor-Disc®, MinElute® (QIAGEN Group); PAXgene® (PreAnalytix GmbH); Roche®, LightCycler® (Roche Group); ABI PRISM®, Applied Biosystems®, StepOnePlus™, ViiA™, ROX™, SYBR® (Life Technologies Corporation); Eppendorf®, Mastercycler® (Eppendorf AG); Stratagene®, Mx3005P®, Mx3000P®, Mx4000® (Agilent Technologies); Bio-Rad®, iCycler®, Chromo4™, CFX96™, DNA Engine Opticon®, CFX384™, iQ™, MyiQ™ (Bio-Rad Laboratories, Inc.); Excel® (Microsoft, Inc.). Registered names, trademarks, etc. used in this document, even when not specifically marked as such, are not to be considered unprotected by law.

HB-0903-003 © 2018 QIAGEN, all rights reserved.

Ordering www.qiagen.com/shop | Technical Support support.qiagen.com | Website www.qiagen.com